

Jan 10, 2026

Cloning with restriction enzymes protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.261ge5krog47/v1>

Julia Heiby¹

¹FLI Leibniz Institute on Aging



Julia C. Heiby

FLI Leibniz Institute on Aging

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.261ge5krog47/v1>

Protocol Citation: Julia Heiby 2026. Cloning with restriction enzymes protocol. protocols.io
<https://dx.doi.org/10.17504/protocols.io.261ge5krog47/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 22, 2024



Last Modified: January 10, 2026

Protocol Integer ID: 106178

Keywords: using restriction enzyme, restriction enzymes protocol, restriction enzyme, gene fragment, gene fragment of interest, enzyme

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

How to quickly clone your gene fragment of interest into a vector, using restriction enzymes.



Materials

- Centifuge 5804R (#5805000010, Eppendorf) or 5810R (#5811000015, Eppendorf)
- Heatblock (#SBH130D3 or #SBH200D3, Stuart)
- 1.5ml tubes (Eppendorf, #0030 120.086)
- Serologicla Pipettes (Greiner, cellstar)
- Milli-Q water system (Merck, Advantage A10)
- E.coli DH5 α (Thermo Fisher, EC0112)
- XhoI (Biolabs, #R01468)
- NotI (Biolabs, #R3189L)
- T4 ligase (Biolabs, #M0202T)
- agarose (Roth, #22675)
- Gel extraction kit (QIAEX®II-QIAGEN, #20021)
- DNA clean up kit (Zymo DNA Clean and Concentrate, #D4006)
- ice
- Tryptone (XXX)
- NaCl (XXX)
- Yeast (XXX)
- Ampicillin (1XX)
- Agar (XXX)
- Mini-Prep kit for DNA isolation (QIAprep®spin, 27106)



Generating Gene Fragment and cutting vector

1

Restriction Enzymes depending on your choice, here for example XhoI and NotI (both Biolabs, R01468 and R3189L , respectively), digesting gene fragments for 3h 37°C, or following the enzyme manufacturer's instruction

+1µl XhoI
+1µl NotI
+2µl cut smart buffer
+ Xµl gene fragment to be cut
Up to 20µl with ddH₂O

Note: If not sure which enzymes to use, check your sequence, for example with SnapGene

DNA clean up

2

Either run the digested fragment on 2% agarose gel (Roth, 22675), cut out bands and use a gel extraction kit (QIAEX®II-QIAGEN, 20021), or, use directly DNA clean up kit by for example Zymo DNA Clean and Concentrate (D4006) and follow manufacturer's instruction

Ligation

3

Ligation with T4 ligase (Biolabs) at 16°C o/n
+ 2µl T4 ligase buffer
+ 1µl T4 ligase
+ Xµl cut gene fragment
+ Xµl cut vector
Up to 20µl with ddH₂O

Note: to calculate the ratio needed between gene fragment and vector, we recommend using the NEB® Calculator, as the amount required will vary depending on the length of gene fragment and vector

Transformation

4

Transformation into competent E.coli DH5α (Thermo Fisher, EC0112)
+ 5µl ligation sample
+ 40µl E.coli DH5α



- incubate for 30min on ice
- heat shock at 42°C for 1min
- 5min on ice
- 1h in LB media 1ml at 37°C
- centrifuge and discard ~900µl of supernatant
- resuspend in leftover media and plate ~50-100µl on LB-plates containing antibiotic (selected resistance of plasmid)
- incubate at 37°C o/n

Amplification

- 5 pick single colonies from plate and transfer to 5ml LB-medium containing antibiotic of choice
incubate at 37°C on shaker o/n

Note:

LB-medium composition: MilliQ H₂O (filled to 1l, and adjusted pH 7.4), Tryptone (10g), NaCl (10g) and Yeast (5g); autoclaved and added for example Ampicillin (100mg/ml stock, add 1ml to 1l LB-medium) when broth is cooled to ~50°C;

Same composition for LB-plates, with addition of agar (15g for 1l) before autoclaving

DNA extraction

- 6 use Mini-Prep kit for DNA isolation following the manufacturers instructions (QIAprep®spin, 27106)

Sequence validation

- 7 Sequence plasmid with appropriate primers and check sequencing results