



May 15, 2024

CODEX FFPE Microtome Sectioning

DOI

<https://dx.doi.org/10.17504/protocols.io.n2bvjn835gk5/v1>

Joanna Bi¹

¹Stanford University



Joanna Bi

Stanford University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n2bvjn835gk5/v1>

Protocol Citation: Joanna Bi 2024. CODEX FFPE Microtome Sectioning. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.n2bvjn835gk5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 15, 2024



Last Modified: May 15, 2024


Protocol Integer ID: 99815

Keywords: codex ffpe microtome sectioning protocol, sectioning ffpe block, ffpe blocks for use, codex

Abstract

Protocol for sectioning FFPE blocks for use with CODEX

- 1 Place the paraffin blocks on wet ice for 5-10 minutes to rehydrate
- 2 Heat up a water bath to 40° C
- 3 Insert a new blade into the microtome
- 4 Prepare Superfrost Plus Microscope slides (Cat. No: 1255015) next to the microtome
- 5 Place a paraffin block on to the microtome and cut into it until the desired part of the sample is reached
- 6 Section the tissue at 4 μm
- 7 Place the sectioned tissue in the water bath until it flattens out
- 8 Place a prepared slide in the water bath, move it towards the tissue, and gently lift it out of the bath
- 8.1 Make sure that the tissue is within these boundaries

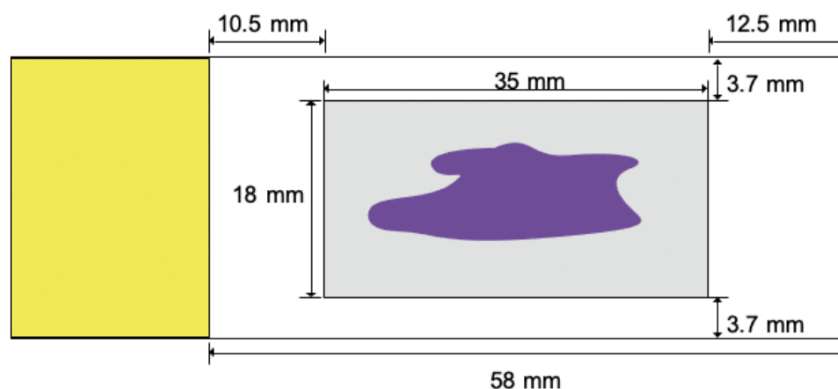



- 9 Change the blade (or move it so an unused portion is being used) between samples and when there is paraffin build up or chipping on the blade
- 10 Let the slides dry facing upwards at an angle
- 11 Store the slides at 4° C for up to 6 months