



Jun 30, 2023

Version 2

Collecting needle and branch samples for terpenoid and expression analysis V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzbryrvx1/v2>

Vidya Vuruputoor¹, karlfetter¹, victoriaburton¹, Jill Wegrzyn¹


¹UConn

PlantCompGenomics



Vidya S Vuruputoor

University of Connecticut

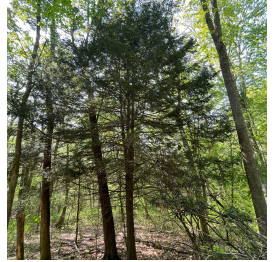

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzbryrvx1/v2>

Protocol Citation: Vidya Vuruputoor, karlfetter , victoriaburton , Jill Wegrzyn 2023. Collecting needle and branch samples for terpenoid and expression analysis. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.rm7vzbryrvx1/v2> Version created by **Vidya S Vuruputoor**



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 25, 2023

Last Modified: June 30, 2023

Protocol Integer ID: 82480

Keywords: branch samples for terpenoid, branch tissue from hemlock conifer, terpenoid analysis, pcg group for sampling needle, hemlock conifer, branch tissue, terpenoid, branch sample, rna expression analysis, sampling needle, pcg group, collecting needle, expression analysis this protocol, expression analysis, sample

Abstract

This protocol is used by the PCG group for sampling needle and branch tissue from hemlock conifers for terpenoid analysis, as well as RNA expression analysis.

Image Attribution

The image was taken by Dr. Karl Fetter



- 1 Open Treesnap (<https://treesnap.org/>), record the accession number on the tree and fill in notes for hemlock woolly adelgid infestation (HWA) levels (scale provided by the app).



3:41

70

<

Hemlock

?

ADD ENTRY

INFORMATION

Images

3 photos added



Species

Eastern hemlock (*Tsuga canadensis*)

⌵

What percentage of the branches you see have hemlock Woolly adelgids?

 SHOW EXAMPLES

☐ 0%

☐ 1-24%

☐ 25-49%

☐ 50-74%

☐ 75-100%

☐ I'm not sure

CANCEL

Canopy health

1 - Healthy

⌵

Tree diameter


5 cm

⌵

Treated

Yes

⌵

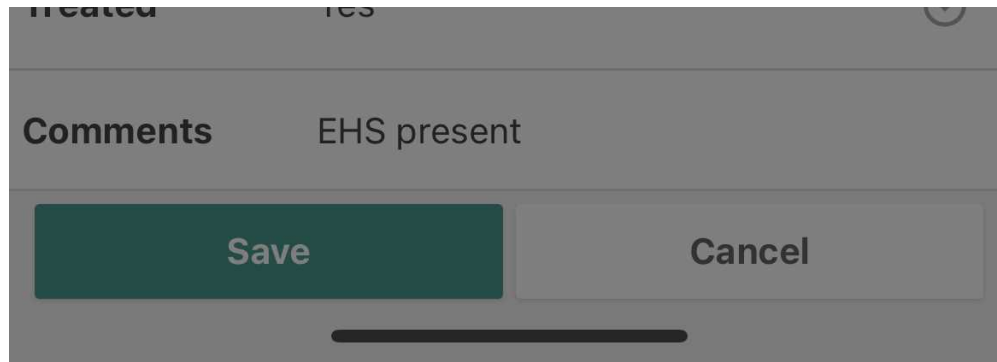

- 1.1 Use the comment field to enter: *EHS present* OR *EHS absent* to record the hemlock scale.

3:40

Hemlock

ADD ENTRY

INFORMATION

Woolly adelgids

1 photo added

Cones

No

Crown classification

Overtopped. This tree's crown is entirely below other trees nearby.

Habitat

Roadside, urban, suburban, or park

Trees nearby

Healthy and large

Canopy health

1 - Healthy

Tree diameter

5 cm

Treated

Yes

Comments

EHS present

Tree Identifier

335-2011E

Advanced Options

**Location**

42.29700, -71.12577

Accuracy 4 meters

[More Options](#)**Save****Cancel**

- 1.2 Take at least 3 photos with TreeSnap. This includes at least one of the trees itself, one photo of the tree tag if that is present, and one photo of the collection vial with the label as well.
- 1.3 Geolocation will be recorded automatically but TreeSnap can also connect with a more precise external GPS device.
- 1.4 Note DBH in appropriate column
- 1.5 Note any other observation you may have- including plant condition (Excellent/Good/Poor), and placement of individual (lower canopy, full sun, etc.)
- 2 If gathering samples for pooled data, choose four random areas around the tree to collect needle and branch samples

Note

For RNA collections- please use gloves, and spray RNase before collections (separate vial)

- 3 Look for the second branch from the shoot tip (stick to the right first) and snip off 5 cm to place in 10 ml cryovial
- 4 Repeat for the other 3 directions (check appropriate column in the sheet provided). Place about 3-4 stems in each 10 ml cryovial.
- 5 Close cryovial, and using the appropriate gloves to handle liquid nitrogen (LN), place cryovial in LN dewar.



Safety information

Use appropriate wear while handling liquid nitrogen

- 5.1 Cap vial tightly to prevent opening in the LN.
- 5.2 Depending on dewer configuration, you can tie tubes with string to a 'halo' to help keep vials from falling to the bottom. Fully submerged tubes can open.

Note

You can hold upto 12 cryovials in a 6L dewer

- 6 Clean trimmers between trees with ethanol (spray bottle).

Note

For collection for RNA- spray clippers with RNase as well

- 7 Use clipboard to manually record notes on a pre-printed spreadsheet with sample names.
- 8 After returning from the field, ensure that notes are consistent with TreeSnap entries. These entries can be amended after collection as needed.