

Apr 30, 2018

Version 1

Colony PCR on *V. natriegens* V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.pukdnuw>

Daniel Bauersachs¹

¹Philipps-Universität Marburg



Daniel Bauersachs

Philipps-Universität Marburg

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.pukdnuw>

Protocol Citation: Daniel Bauersachs 2018. Colony PCR on *V. natriegens*. protocols.io

<https://dx.doi.org/10.17504/protocols.io.pukdnuw>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 30, 2018



Last Modified: April 30, 2018

Protocol Integer ID: 11884

Keywords: colony pcr, pcr



1

5 µl	reaction buffer
0,5 µl	dNTPs
0,5 µl	primer foward
0,5 µl	primer reverse
0,1 µl	one-taq polymerase
18,4 µl	A.dest.
= 25 µl	

Now you have to pick one single colony with a pipette tip and put it into your reaction mix. After ~15 seconds you can take the tip out and start the PCR.