

Sep 23, 2019

Colony PCR (Protocol for Thermo Scientific™ Phire™ Hot Start II DNA Polymerase)

DOI

dx.doi.org/10.17504/protocols.io.7kkhkuw

Alba Balletbó¹

¹Wageningen University

iGEM Wageningen 2019



Alba Balletbó

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.7kkhkuw

Protocol Citation: Alba Balletbó 2019. Colony PCR (Protocol for Thermo Scientific™ Phire™ Hot Start II DNA Polymerase). [protocols.io](https://dx.doi.org/10.17504/protocols.io.7kkhkuw) <https://dx.doi.org/10.17504/protocols.io.7kkhkuw>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 23, 2019

Last Modified: September 23, 2019

Protocol Integer ID: 28012



Abstract

Thermo Scientific™ Phire™ Hot Start II DNA Polymerase is a novel DNA polymerase designed for use in all routine and high throughput PCR applications. A special DNA binding domain linked to the Phire Hot Start II DNA Polymerase enhances the processivity of the polymerase, enabling short extension times and improved yields. The polymerase is also capable of amplifying long DNA fragments, such as the 7.5 kb genomic DNA used in Thermo Scientific quality control assays. Phire Hot Start II DNA Polymerase provides 2-fold fidelity compared to Taq DNA polymerase. The hot start modification of the polymerase is based on the Affibody™ inactivation method^{1,2}. It inhibits DNA polymerase activity at ambient temperatures, thus preventing amplification of non-specific products. At polymerization temperatures the Affibody molecule is released, rendering the polymerase fully active. Phire Hot Start II DNA Polymerase generates blunt ends in the amplification products. It does not possess the 5'→3' exonuclease activity needed for hydrolysis experiments.

Guidelines

The annealing rules are different from many common DNA polymerases (such as Taq DNA polymerases).

Use 10–15 s/kb for extension.


Phire Hot Start II DNA Polymerase produces blunt end DNA products.

Materials

MATERIALS

 Phire Hot Start II PCR Master Mix **Thermo Fisher Catalog #F125S**

Safety warnings

 Wear laboratory coat and glasses during the whole procedure.

- 1 Add all components in a 250 μL tube making up to a 20 or 50 μL reaction.

'Pro-tip': Carefully mix and centrifuge all tubes before opening to ensure homogeneity and improve recovery. When using Phire Hot Start II PCR Master Mix, it is not necessary to perform the PCR setup on ice.

'Note': Due to the unique nature of Phire Hot Start II DNA Polymerase, optimal reaction conditions may differ from standard enzyme protocols.

'Pro-tip': Phire Hot Start II DNA Polymerase tends to work better at elevated denaturation and annealing temperatures due to higher salt concentrations in its buffer.

Add items in this order:

Com pone nts	20 μL reacti on	50 μL reacti on	Final conc entra tion
H ₂ O	add to 20 μL	add to 50 μL	
2X Phire Hot Start II PCR Mast er Mix	10 μL	25 μL	1X
Forw ard prime r *	X μL	X μL	0.5 μM
Rever se prime r *	X μL	X μL	0.5 μM
Temp late DNA	X μL	X μL	
DMS O**, optio nal	0.6 μL	1.5 μL	3%

* The recommendation for final primer concentration is 0.5 μM , but it can be varied in a range of 0.2–1.0 μM , if needed.

** Addition of DMSO is recommended for GC-rich amplicons. DMSO is not recommended for amplicons with very low GC %.

- 2 Gently mix the PCR reactions and transfer the tubes to a thermocycler. Thermocycling conditions for a routine PCR:

	Step	Temp eratu re	Time	Cycle s
	Initial Denat uratio n	98°C	30 s	1
	Denat uratio n	98°C	5 s	25- 35
	Anne aling*	X°C	5 s	
	Exten sion* *	72°C	10-15 s/kb	
	Final Exten sion	72°C	1 min	1
	Hold	4°C		

* The optimal annealing temperature for Phire Hot Start II DNA Polymerase may be significantly different than annealing temperature with other DNA polymerases. Always use the T_m calculator and instructions from the supplier to determine the T_m values of the primers and optimal annealing temperature. As a basic rule, for primers > 20 nt, anneal for 5 seconds at a T_m +3 °C of the lower T_m primer. For primers ≤ 20 nt, use an annealing temperature equal to the T_m of the lower T_m primer. If necessary, use a temperature gradient to find the optimal annealing temperature for each template-primer pair combination.

** The extension should be performed at 72 °C. Extension time of 10 seconds per 1 kb is recommended for most templates. However, higher yields may be obtained using extension time of 15 s/kb with challenging primer template pairs.