

Sep 09, 2020

## Colorimetric RT-LAMP Research group SARS-COV-2 detection protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.bk59ky96>

Teodor Emilov Mentov<sup>1</sup>

<sup>1</sup>HUBU

XPRIZE Rapid Covid Tes...



vnogi

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



**DOI:** <https://dx.doi.org/10.17504/protocols.io.bk59ky96>

**Protocol Citation:** Teodor Emilov Mentov 2020. Colorimetric RT-LAMP Research group SARS-COV-2 detection protocol. protocols.io <https://dx.doi.org/10.17504/protocols.io.bk59ky96>

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

We use this protocol and it's working

**Created:** September 09, 2020

**Last Modified:** September 09, 2020

**Protocol Integer ID:** 41889

**Keywords:** lamp research group sar, sar, cov, detection

## vortex swab sample

- 1 Vortex swab sample in Viral transport medium
- 2 Take 200µl and transfer to 1,5ml microtube
- 3 Add 500ul of ethanol reconstituted lysis buffer incubate 10 minutes. Vortex or shake after
- 4 Add solubilization buffer, and wait 2 minutes, shake or vortex after
- 5 Add 200µl 98% Ethanol shake and use a an intense and brief centrifuge. 5-10 sec to 30 000g. This will avoid clogging in the filter cartrige afterwards.
- 6 Transfer the supernatant to the Kurabo catrige.
- 7 Wash with wash buffer and filter 3 times with Kurabo-Mini480.
- 8 add 50-100 (better 50)µl of elution buffer or pcr grade water.
- 9 filter and you have extracted the viral RNA in the elution tubes. You may store it in -80C or use it directly in the isothermic amplification
- 10 Isothermal amplification: Make a mix of 5µl of eluate sample and 5µl of pcr grade H2O and 2,5 µl of 10x primer mix in a 1,5ml transparent microtube. Standart primer concentrations of 1·6 µM FIP/BIP, 0·2 µM F3/B3, 0·4 µM FL/BL work just fine.
- 11 add 12.5 µl of colorimetric RT-LAMP master mix. The NEB one or your own. Mix well with the pipette tip.
- 12 add 50µl of mineral oil to seal the reaction. This is essencial. Unsealed reactions are very difficult to detect.



- 13 Make a photo with your mobile phone. to register base colour, should be red-pink.
- 14 incubate 30 minutes at 63C°. Take a photo afterwards. You dshold see the positive reactions as yellow. You can register it measuring the green channel in the RG colourspace (a normalized RGB colourspace). Higher than 83 is positive, loyer or equal is negative. To do this take the median values of the RGB measure with ImageJ, divide R,G,B to R+G+B. That is RG colorspace. You only need the green channel.