



Oct 10, 2018

Confocal Microscopy imaging for Opaline Silica Single Cell Skeletons (Polycystines Radiolaria)

DOI

<https://dx.doi.org/10.17504/protocols.io.t5qeq5w>

Miguel Mendez Sandin¹

¹Station Biologique de Roscoff

Ecology of Marine Plankton (ECOMAP) team - Roscoff
Tech. support email: phytopk@sb-roscoff.fr



Miguel Mendez Sandin

Institute of Evolutionary Biology (CSIC-UPF)

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.t5qeq5w>

Protocol Citation: Miguel Mendez Sandin 2018. Confocal Microscopy imaging for Opaline Silica Single Cell Skeletons (Polycystines Radiolaria). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.t5qeq5w>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: September 30, 2018

Last Modified: October 10, 2018

Protocol Integer ID: 16272

Keywords: opaline silica single cell skeleton, confocal microscopy imaging, microscopy, polycystines radiolaria, single cell, radiolaria

Recovering skeletons

- 1 After DNA extraction, recover skeleton from the eluted pellet under binoculars or inverted microscope.

Note: During DNA extraction: dilute waste from the extraction procedure (i.e. pellet debris, containing the skeleton) in milliQ water and store skeletons at -20°C.

Rinsing

- 2 Rinse skeleton several times in milli-Q water to decrease the concentration of SDS and other lysis and DNA precipitation reagents.

Cleaning

- 3 Transfer skeleton into 1.5 ml Eppendorf tubes containing 50 µl of hydrogen peroxide (H₂O₂).
- 4 Heat at 70°C for 10 min to remove residual organic matter.

Diluting

- 5 Add 1 ml of milli-Q water.

Rinsing

- 6 Handpick skeleton under binoculars or inverted microscope and repeat several rinsing steps.

Staining

- 7 Transferred into 1.5 ml Eppendorf tubes containing 25µl of fluorescent staining mix (Colin et al., 2017).



Colin, S., Coelho, L.P., Sunagawa, S., Bowler, C., Karsenti, E., Bork, P., Pepperkok, R., de Vargas, C., 2017. Quantitative 3D-imaging for cell biology and ecology of environmental microbial eukaryotes. *Elife* 6, 1–15. doi:10.7554/eLife.26066

- 8 Let the tubes for two hours for skeleton staining at room temperature and dark conditions.

Diluting

- 9 Add 1.5 ml of milli-Q water to dilute the dye.

Rinsing

- 10 Handpick the skeleton and repeat 4-6 rinsing steps into milli-Q water to properly clean the sample.

Preparing for imaging

- 11 Transfer skeleton into a micro well plate.
- 12 Dry at 70 °C for 2 hours.

Imaging

- 13 Skeleton ready to be imaged