



Jun 21, 2023

Version 1

🌐 Converting ssDNA oligos to dsDNA with T4 DNA polymerase V.1

DOI

dx.doi.org/10.17504/protocols.io.261ged6xyv47/v1

Alex N Nguyen Ba¹

¹University of Toronto



Alex N Nguyen Ba

University of Toronto

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.261ged6xyv47/v1>

Protocol Citation: Alex N Nguyen Ba 2023. Converting ssDNA oligos to dsDNA with T4 DNA polymerase. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.261ged6xyv47/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 21, 2023



Last Modified: April 18, 2024

Protocol Integer ID: 83806

Keywords: ssdna to dsdna oligo, t4 dna polymerase this protocol, t4 dna polymerase, dsdna oligo, converting ssdna, ssdna, fidelity polymerase, polymerase, dna, complementary oligo, order of the extension temperature, high temperature of typical thermocycling

Abstract

This protocol allows one to convert ssDNA to dsDNA oligos. In principle, one can buy two complementary oligos and anneal them. However, there are a few cases where randomized bases are desired, and therefore complementary oligos can not be ordered. The protocol works similarly to PCR, except that the reaction is not performed at the high temperature of typical thermocycling, because the oligos are short and their melting temperature can be in the order of the extension temperature of high-fidelity polymerases.

Protocol materials

☒ NEBuffer 2.1 - 5.0 ml **New England Biolabs Catalog #B7202S**

☒ dNTPs

☒ T4 DNA Polymerase - 750 units **New England Biolabs Catalog #M0203L**

☒ EDTA (0.5 M), pH 8.0 **Life Technologies Catalog #AM9260G**

☒ Monarch PCR and DNA Cleanup Kit - 250 preps **New England Biolabs Catalog #T1030L**


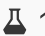


Troubleshooting



Primer design



- 1 Design your target single-stranded DNA oligo such that the 3' end contains the reverse complement of an extending oligonucleotide. For example, appending "GTCATAGCTGTTTCCTG" to the end of your oligo will allow an oligonucleotide matching the M13 Reverse (-27) primer to extend it (5'-CAGGAAACAGCTATGAC-3').
- 2 Resuspend your oligonucleotides in 1/10th TE (10mM Tris, 0.1mM EDTA pH 8) to a final concentration of 100uM to form your stock oligonucleotide solutions.
- 3 Make working oligonucleotide solutions by diluting the stock oligonucleotides to a final concentration of 10uM.

Oligo annealing

- 4 Mix in a PCR tube  5 μL of a [M] 10 micromolar (μM) stock of the target oligo with  10 μL of a [M] 10 micromolar (μM) stock of the extending oligo. Add 27ul of molecular biology grade water.
- 5 Add  5 μL of  NEBuffer 2.1 - 5.0 ml **New England Biolabs Catalog #B7202S** . Mix well by vortexing.
- 6 Place the tube in a thermocycler and run the following protocol:
 1. 95 degrees for 10 seconds
 2. Decrease by 1 degrees
 3. Repeat step 2 every 10 seconds, for 90 cycles.
 4. Hold at 4 degrees.

15m

Oligo extension



- 7 Add  2.5 μL  dNTPs ([M] 10 millimolar (mM) stock) and mix well by vortexing.
- 8 Go to the thermocycler and prepare the reaction cycle. When ready, start the protocol and pause when the block reaches 0 degrees.
 1. 5 minutes at 0 degrees C.
 2. 5 minutes at 22 degrees C.

40m



3. 30 minutes at 37 degrees C.



4. Hold at 0 degrees.

- 9 Add  0.5 μ L of  T4 DNA Polymerase - 750 units **New England Biolabs Catalog #M0203L** and mix well.

- 10 Place the tube in the thermocycler and unpause the run.


Oligonucleotide cleanup

10m

- 11 As soon as the reaction is completed, add  EDTA (0.5 M), pH 8.0 **Life Technologies Catalog #AM9260G** to  10 millimolar (mM) .

- 12 Use a PCR purification kit to purify the oligo. Ensure the kit is able to purify short oligonucleotides if needed (a kit such as

10m

 Monarch PCR and DNA Cleanup Kit - 250 preps **New England Biolabs Catalog #T1030L**

can be suitable).