

Aug 24, 2020

Copy of ELISA for quantification of human C5 in serum or plasma.

DOI

dx.doi.org/10.17504/protocols.io.bj76krre

Angel A Justiz-Vaillant¹, Belkis Ferrer-Cosme²

¹University of the West Indies St. Augustine;

²"Saturnino Lora Torres' Provincial Teaching Clinical Surgical Hospital. Cuba



Angel A Justiz-Vaillant

University of the West Indies St. Augustine

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bj76krre

Protocol Citation: Angel A Justiz-Vaillant, Belkis Ferrer-Cosme 2020. Copy of ELISA for quantification of human C5 in serum or plasma.. [protocols.io https://dx.doi.org/10.17504/protocols.io.bj76krre](https://dx.doi.org/10.17504/protocols.io.bj76krre)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: August 24, 2020

Last Modified: August 24, 2020

Protocol Integer ID: 40926

- 1 An anti-human C5 coating antibody is adsorbed onto the microwells by incubation overnight at 4°C with carbonate-bicarbonate buffer.
- 2 Add 50 µl of human serum or plasma. Human C5 present in the serum or plasma binds to antibodies adsorbed into the microwells.
- 3 The microplate is blocked with 3% non-fat milk-PBS buffer and later wash to remove unbound proteins.
- 4 Fifty (50) µl of biotin-conjugated anti-C5 antibody is added. The optimal dilution must be investigated.
- 5 The microplate is rewashed with PBS-Tween 20 buffer, pH 7.4.
- 6 One hundred µl of streptavidin-HRP conjugate is added and it binds to the biotin-conjugated anti-C5 antibody. The optimal dilution of this conjugate must be investigated.
- 7 The plate is washed following incubation to remove the unbound Streptavidin-HRP.
- 8 Add 100 µl of 3,3',5,5'- tetramethylbenzidine (TMB; Sigma-Aldrich) into each well.
- 9 Incubate the microwells in the dark for 15 min.
- 10 A colored product is formed in proportion to the quantity of C5 present in the sample or standard.
- 11 The reaction is terminated by addition of 100 µl 3M H₂SO₄ and the absorbance is measured at 450 nm.
- 12 A standard curve is made from 7 human C5 standard dilutions and the human C5 sample concentration is determined.