

Nov 28, 2019

## Copy of Protocol for the production of crude alcoholic extracts from native plants

DOI

[dx.doi.org/10.17504/protocols.io.9vch62w](https://dx.doi.org/10.17504/protocols.io.9vch62w)



Marissol M Leite<sup>1</sup>, Hadda Tercya Lima Silva<sup>2</sup>, Zanderluce Gomes Luis<sup>2</sup>, Aline Coutinho Cavalcante<sup>2</sup>, Caio Maximino<sup>2</sup>, Diogenes Silva<sup>2</sup>

<sup>1</sup>Universidade federal do sul e sudeste do Pará; <sup>2</sup>Universidade Federal do Sul e Sudeste do Pará

Medicinal Plants Southe...



Diogenes Silva

Universidade Federal do Sul e Sudeste do Pará, Universidade ...

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.9vch62w>

**Protocol Citation:** Marissol M Leite, Hadda Tercya Lima Silva, Zanderluce Gomes Luis, Aline Coutinho Cavalcante, Caio Maximino, Diogenes Silva 2019. Copy of Protocol for the production of crude alcoholic extracts from native plants. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.9vch62w>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** November 28, 2019

**Last Modified:** November 28, 2019

**Protocol Integer ID:** 30340

**Keywords:** amazonian plants, crude extract, fish anesthetize, crude alcoholic extracts from plant organ, crude alcoholic extracts from native plant, obtaining crude alcoholic extract, production of crude alcoholic extract, crude alcoholic extract, native plants this protocol, used plant, piper callosum, native plant, spilanthes acmella, zebrafish danio rerio, plant, plant organ, such leaf, joão brandinho

## Abstract

This protocol describes the method to obtaining crude alcoholic extracts from plant organs, such leaves and inflorescences. The used plants were the popularly called "João Brandinho" *Piper Callosum* and "Jambu" *Spilanthes acmella*. The crude extracts were used in our laboratory to anesthetize the native silver tetra fish *Ctenobrycon sp.* and Zebrafish *Danio rerio*, by immersion method.

## Materials

### MATERIALS

☒ Distilled Water

☒ 70% alcohol Merck MilliporeSigma (Sigma-Aldrich) Catalog #793213

☒ Paper towels

☒ Beaker

☒ Fluted Qualitative Filter Paper Circles Thermo Fisher Catalog #0979014G

☒ Wyllie Micro - te 650 Knife Mill

☒ Analytical Balance

☒ Amber glass

☒ Tray

☒ Glass funnel

☒ crucible

## Troubleshooting

- 1 Separate the leaves and inflorescences from plants with the help of scissors.



- 2 Rinse the material under running water and then with distilled water to remove debris.



- 3 Dry the material with paper towels to remove the excess of water.



- 4 Grind the leaves and flowers with a knife mill or blender and / or just cut with the help of scissors, in order to allow a better extraction of secondary compounds.





- 5 Weight the vegetable mass into an analytical balance and dilute in 70% alcohol.
- 6 After quantifying the vegetable mass, dilute and mash the crushed and / or cut material into 70% alcohol in the desired ratio, such as 1:1 (1 gram of vegetal material to 1 milliliter of 70% alcohol) until the thoroughly solution mix.



- 7 After maceration and dilution in 70% alcohol filter the obtained solution with the aid of filter paper in a beaker.



- 8 Identify the alcoholic crude extract and store in an amber glass in the refrigerator