

Jul 16, 2024

Coronal cryosectioning of mouse brains

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g713w9gwz/v1>

Lauren Faget¹, Thomas Hnasko²

¹UCSD; ²University of California, San Diego



Lauren C. Faget

Boise State University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g713w9gwz/v1>

Protocol Citation: Lauren Faget, Thomas Hnasko 2024. Coronal cryosectioning of mouse brains. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.q26g713w9gwz/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 16, 2024



Last Modified: August 13, 2024

Protocol Integer ID: 103431

Keywords: coronal cryosectioning of mouse brains cryostat, mouse brains cryostat, coronal cryosectioning, mouse brain tissue, mouse brain, brain

Disclaimer

The [protocols.io](https://www.protocols.io) team notes that research involving animals and humans must be conducted according to internationally-accepted standards and should always have prior approval from an Institutional Ethics Committee or Board.

Abstract

Cryostat coronal cryosectioning of mouse brain tissue

Before start

PFA-perfused, 30% sucrose-cryoprotected, and flash-frozen mouse brains are stored at -80C until cryosectioning

Procedure

25m

- 1 Retrieve frozen mouse brain stored at  -80 °C .
- 2 Place brain in cryostat. Set chamber temperature to  -16 °C to  -20 °C . Set brain holder temperature to  -20 °C . Let brain equilibrate to temperature for at least  00:20:00 before sectioning. 20m
- 3 Insert chuck and blade into cryostat.
- 4 Apply a drop of OCT on the chuck, and affix brain vertically (olfactory bulbs up).
- 5 Apply another layer of OCT at the base of the brain, turning the brain-holder with the other hand. Cover the cerebellum and the junction between the cerebellum and cortex with OCT. Allow OCT to freeze solid (~  00:05:00). 5m
- 6 Cut 16 to 30 micron sections. Adjust angle of medio-lateral and ventro-dorsal axis so that sections are free of cutting artifacts, symmetrical, and well-aligned to reference atlas.
- 6.1 Use of anti-roll plate is optional and can help reduce rolling of sections, but must be carefully adjusted to avoid cutting artifacts.
- 6.2 For immunohistochemistry we typically cut 30-micron sections for best results. For RNAscope we typically cut 16 to 20-micron sections.
- 7 Collect free-floating sections in 48-well plate.
- 7.1 Label your plate lid and base with brain/experiment details using lab marker and/or lab tape.
- 7.2 Pre-fill wells with  1 mL 1X PBS containing 0.01% sodium azide (100mg/L).



- 7.3 Can use syringe with blunt needle to collect sections after sectioning. Dip needle into PBS and collect section by gently touching it on the side of the section.
- 7.4 Collect 5 sections per well. Or for experiments that require finer resolution of coordinate location can collect sections serially, adding 1 section per well until the end of the plate and start back at well #A1.
- 8 Use parafilm to seal the lid to the plate (air-tight) and refrigerate at  4 °C until further processing.