



Mar 21, 2024

Correlated light-electron microscopy (CLEM) and cryo-focused ion beam (FIB) milling

DOI

<https://dx.doi.org/10.17504/protocols.io.dm6gpzexjlzp/v1>

Dorothy zhao¹

¹Max planck institute of biochemistry



Dorothy zhao

Max planck institute of biochemistry

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.dm6gpzexjlzp/v1>

Protocol Citation: Dorothy zhao 2024. Correlated light-electron microscopy (CLEM) and cryo-focused ion beam (FIB) milling. [protocols.io https://dx.doi.org/10.17504/protocols.io.dm6gpzexjlzp/v1](https://dx.doi.org/10.17504/protocols.io.dm6gpzexjlzp/v1)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: March 21, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 97061

Keywords: ASAPCRN, interactions of autophagosome, autophagosome, aggregation of protein, amyloid fibril, interactions between the autophagy receptor p62, autophagy receptor p62, autophagy, trapping of autophagic structure, autophagic structure, limited efficiency of autophagy, organized polyq fibril, fibrillar disease aggregate, protein of hd, polyq aggregates in cultured cell, polyq fibril, disease protein of hd, electron microscopy, amyloid, neurodegenerative disease, protein, expanded polyglutamine, disease protein, amorphous aggregate phase, like fibrils by liquid, cytopathologic hallmark, inherited neurodegenerative disease, like fibril, aggrephagy, cell, efficiency of aggrephagy

Funders Acknowledgements:

ASAP

Grant ID: CRN

Abstract

Aggregation of proteins containing expanded polyglutamine (polyQ) repeats is the cytopathologic hallmark of a group of dominantly inherited neurodegenerative diseases, including Huntington's disease (HD). Huntingtin (Htt), the disease protein of HD, forms amyloid-like fibrils by liquid-to-solid phase transition. Macroautophagy has been proposed to clear polyQ aggregates, but the efficiency of aggrephagy is limited. Here, we used cryo-electron tomography to visualize the interactions of autophagosomes with polyQ aggregates in cultured cells *in situ*. We found that an amorphous aggregate phase exists next to the radially organized polyQ fibrils. Autophagosomes preferentially engulfed this amorphous material, mediated by interactions between the autophagy receptor p62/SQSTM1 and the non-fibrillar aggregate surface. In contrast, amyloid fibrils excluded p62 and evaded clearance, resulting in trapping of autophagic structures. These results suggest that the limited efficiency of autophagy in clearing polyQ aggregates is due to the inability of autophagosomes to interact productively with the non-deformable, fibrillar disease aggregates.

CLEM

- 1 Vitrified sample on autogrids were loaded onto a cryo-confocal LM set up (Leica SP8) equipped with a 50X/0.9 NA objective (Leica Objective), metal halide light source (EL6000), air-cooled detector (DFC900GT), a cryo-stage (-195 °C), and two HyD detectors.
- 2 The sample was kept in liquid nitrogen vapor, following a similar workflow. Avoid keeping sample in instrument for extended time due to high ice contamination. Best to finish a grid within 30 minutes.
- 3 Cryo-confocal z-stacks (step size 500 nm, x-y pixel size 85 nm) were taken with pin hole = 1, and a 9 µm depth with the LAS X Navigator software, using 488 and 552 nm laser excitation for GFP and RFP tagged proteins, respectively, also picking up signals from the auto-fluorescent Dynabeads. Depending on target size, the depth and increment of the scan could be changed for optimization.
- 4 To improve signal clarity, image stacks were de-convoluted and restored with Huygens Essential software (Scientific Volume Imaging) to remove noise.
- 5 The stack was then imported into the 3Dcorrelation software (Arnold et al., 2016) and re-sliced into cubic voxels.

FIB milling

- 6 For preparing the lamella (150-250 nm), autogrids were mounted into a Quanta dual-beam 3D FIB/scanning **electron microscope** (Thermo) equipped with a transfer shuttle system (PP3000T, Quorum) at < -180 °C throughout milling.
- 7 To protect the milling front of the lamella, gaseous organometallic platinum was sprayed onto the sample on the cryo-stage using a gas injection system. The duration of the gas spray varies between instruments.
- 8 To target the cell for milling, the grid square was correlated with the cryo-confocal fluorescence z-stack using the 3Dcorrelation software. The target was imaged and correlated iteratively with the z-stack throughout milling for accuracy. One can check for accuracy by correlating the SEM image with the fluorescence image if the fluorescent target is large.



- 9 The 12-15 μm wide lamellas were generated using a Gallium FIB at 30 kV with a 20° stage angle in three consecutive steps. The more distant region ($>2\text{ }\mu\text{m}$) above and below the target was rough milled with a higher current of 500 pA, followed by fine milling to a $\sim 800\text{ nm}$ lamella using a current of 100 pA.
- 10 A final polishing of the lamella to thickness of 150-250 nm was carried out with a 30-50 pA current. Lamella final thickness was estimated with SEM at 3 keV, for a lack of overcharging.