



Jun 26, 2023

Cortical, Striatal and Dopaminergic Neurons Cultures Protocol

DOI

dx.doi.org/10.17504/protocols.io.rm7vzxn48gx1/v1

Giulia Tombesi¹

¹Northwestern University, Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD 20815



Giulia Tombesi

Northwestern University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzxn48gx1/v1>

Protocol Citation: Giulia Tombesi 2023. Cortical, Striatal and Dopaminergic Neurons Cultures Protocol. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.rm7vzxn48gx1/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: June 26, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 84050

Keywords: ASAPCRN, co-cultures, primary neurons, striatal-cortical neuronal cultures, dopaminergic-cortical neuronal cultures, primary neuronal cultures from postnatal mice, dopaminergic neurons cultures protocol this protocol, dopaminergic neurons cultures protocol, pure cortical neurons culture, primary neuronal culture, cortical neuron, dopaminergic neuron, postnatal mice, neuron, striatal, postnatal, mice

Funders Acknowledgements:

Aligning Science Across Parkinson's [ASAP-020600] through the Michael J. Fox Foundation for Parkinson's Research (MJFF)
Grant ID: ASAP-020600

Abstract

This protocol describes how to obtain primary neuronal cultures from postnatal mice (P0-P1). It can be used to obtain pure cortical neurons cultures or stratal-cortical neurons and dopaminergic-cortical neurons co-cultures.

Materials

Papain Dissociation System from worthington biochemical corporation

Troubleshooting



Before start

Prepare the plates:

- Coat plates with poly-L-Lysine Overnight at room temperature or at least 02:00:00 at 37°C.
- Wash thoroughly with distilled water three times. Let it dry under the hood in the dark.

Prepare the following solutions:

- **Dissection medium:** 10mM HEPES pH7.3 , 0.1% D-glucose, 0.11mg/ml Sodium pyruvate in HBSS.
- **DNase solution:** one vial of DNase in 500 μ L of BME.
- **Papain solution:** one vial of Papain in 5 mL of BME plus 250 μ L of DNase solution.
- **Trypsin inhibitor solution:** 15.5mg/ml in BME. Filter the solution.
- **Stop solution:** 250 μ L of DNase solution, 600 μ L of Trypsin inhibitor solution in 5.4 mL of BME.
- **10/10 solution:** Trypsin inhibitor 10ug/ml and BSA 10ug/ml in BME. Filter the solution.
- **Plating medium:** 1 mL B27 supplement, 125 μ L L-Glutamine, 500 μ L Pen/Strep, 500 μ L N2 supplement in 50 mL Neurobasal.



Dissection

50m

- 1 Dissect the cortices, the striatum and the substantia nigra (SN) in cold dissection medium, carefully removing the meninges

Cells Dissociation

1h 5m

- 2 Snip the tissues into smaller pieces and place them in Papain solution. Use around  500 μL of Papain solution for 5 SN, around  1 mL of Papain solution for 5 striata and around  5 mL of Papain solution for 5-6 cortices.

- 3 Triturate 10 times with a 5ml-pipette

- 4 Incubate at 37°C for  00:40:00 and mix by inversion every  00:10:00

50m

Cells Dissociation

1h 5m

- 5 Triturate 10 times again with a 5ml-pipette

Cells Dissociation

1h 5m

- 6 Spin in swing bucket rotor at 1000rpm for  00:05:00

5m

- 7 Remove and discard the supernatant

- 8 Tap the tubes to move the pellet and add STOP solution. Add  1 mL for 5 SN,  1 mL for 5 striata and around  3 mL for 3 cortices.

- 9 Triturate 3 times with a 5ml-pipette



- 10 Incubate at room temperature for  00:10:00 10m
- 11 Take the supernatant and add to 10/10 solution drop-by-drop (use a volume of 10/10 solution that is equal to the supernatant volume)
- 12 Spin at 800rpm for  00:10:00 10m

Neuron Seeding

- 13 Discard the supernatant and resuspend in Neurobasal complete medium with AraC. Use  1 mL of medium to resuspend 5 striata and 3 cortices.
- 14 Count cortical and striatal cells with trypan blue
- 15 Plate 500000 cells/well in 24wells-plate.
For striatal-cortical co-culture plate a ratio of 1:2 striatum:cortex.
For SN-cortical co-culture plate one well for each SN and around 500000 cortical neurons/well and also add GDNF to the medium.
- 16 Replace half medium about every 7 days by collecting conditioned medium from cells and mixing with fresh complete medium in a ratio of 1:1.