

Apr 20, 2022

Counting kidneys

DOI

<https://dx.doi.org/10.17504/protocols.io.6qpvr6k4pvmk/v1>



Gabriel J. Barrero¹, Abraham Palmer¹, Oksana Polesskaya¹

¹UCSD

CGORD



Gabriel J Barrero

Johns Hopkins University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6qpvr6k4pvmk/v1>

Protocol Citation: Gabriel J. Barrero, Abraham Palmer, Oksana Polesskaya 2022. Counting kidneys. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.6qpvr6k4pvmk/v1>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 11, 2022

Last Modified: April 20, 2022


Protocol Integer ID: 60610

Keywords: counting kidney, kidneys in rat, rat


Abstract

Dissecting and counting kidneys in rats.


- 1 Rats are euthanized as per IACUC guidelines
Spray fur with ethanol. Pick up skin above the sternum, and make a cut through skin and muscle layers.
Don't cut the chest cavity (don't go above the diaphragm).
- 2 Cut skin and muscles from the sternum down to the lower belly. Mostly you will see liver and colon.

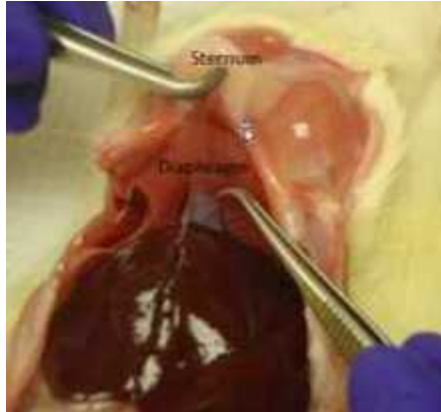

- 3 Lift liver lobes and push aside fat tissue and intestines. You will see dark-colored organs – kidney and spleen.

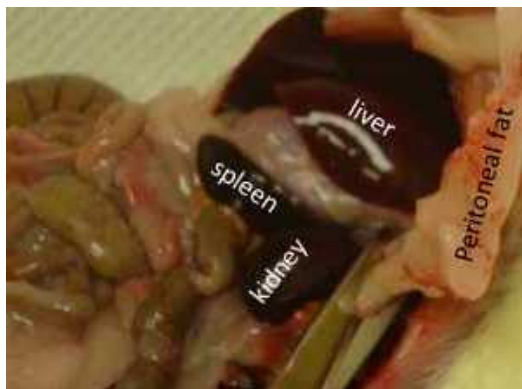

Gross dissection

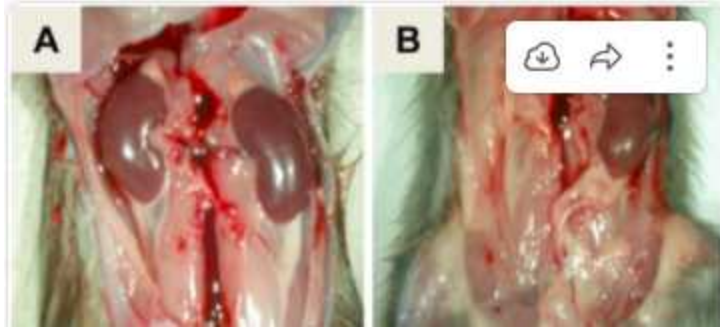

Representative renal anomalies observed in rats. A. Grossly normal renal development. B. Unilateral renal agenesis (URA) PloS one. 10. e0118147. 10.1371/journal.pone.0118147

- 4 Count the number of observed kidneys and make a record in the file.