Aug 27, 2019

## CPEC Protocol



In 1 collection

DOI

dx.doi.org/10.17504/protocols.io.6t7hern

N.J. Hillson<sup>1,2</sup> <sup>1</sup>JBEI; <sup>2</sup>LBNL

**Michael J Fero** 





DOI: dx.doi.org/10.17504/protocols.io.6t7hern

Protocol Citation: N.J. Hillson 2019. CPEC Protocol. protocols.io https://dx.doi.org/10.17504/protocols.io.6t7hern

## **Manuscript citation:**

J. Quan and J. Tian, "Circular polymerase extension cloning of complex gene libraries and pathways," PloS one, vol. 4, no. 7, p. 6441, 2009.

License: This is an open access protocol distributed under the terms of the Creative Commons Attribution License, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working We use these protocols in our group and they work.

Created: August 26, 2019

Last Modified: August 27, 2019

Protocol Integer ID: 27231

- 1 Measure the DNA concentration (ng/ml) of each assembly piece.
- 2 Add 100 ng of the linearized vector backbone and equimolar amounts of the other assembly pieces to a 25 ml total volume assembly reaction mixture as follows:

linearized vector backbone (100 ng)

- + each additional assembly piece (to equimolar with backbone)
- + 5 ml 5X HF Phusion Reaction Buffer
- + 1 ml 10 mM dNTPs
- + 0.75 ml DMSO
- + 0.5 ml 2U/ml Phusion Polymerase

```
+ _____dH<sub>2</sub>O to
```

25 ml

3 Perform the assembly reaction in a thermocycler as follows:

```
30 sec @ 98 C 1 cycle
10 sec @ 98 C }
30 sec @ 55 C } 1 to 15 cycle(s)**
length* (kb) x 15 sec @ 72 C }
10 min @ 72 C 1 cycle
```

\*The total length of the assembled product (in kb) \*\*The number of repeated cycles should exceed the number of assembly pieces

Transform 5 ml of the assembly reaction into 100 ml of competent E. coli and/or run a diagnostic agarose gel to check for successful assembly.