

Feb 21, 2025

CRISPR/Cas9 gene editing in hiPSC

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlk9xxdv5r/v1>

Laura Mahoney¹

¹UCL / The Francis Crick Institute



Laura Mahoney

UCL / The Francis Crick Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlk9xxdv5r/v1>

Protocol Citation: Laura Mahoney 2025. CRISPR/Cas9 gene editing in hiPSC. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.j8nlk9xxdv5r/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 21, 2025



Last Modified: February 21, 2025

Protocol Integer ID: 123201

Keywords: cas9 gene editing in hipsc, cas9 gene editing, cas9 in humain inducible pluripotent cell, point mutation via crisp, crispr, hipsc, cas9, humain inducible pluripotent cell, point mutation, editing

Abstract

This protocol is adapted from Skarnes et al., 2019 and describes the steps to follow for knocking-in point mutation via CRISP/Cas9 in humain inducible pluripotent cells (hiPSC)

A. Pre-nucleofection

- 1 Resuspend the sgRNA overnight at 4C in TE buffer at a concentration of **1M 4 µg/µL**
- 2 Resuspend the Alt-R HDR modified ssODN overnight in D-PBS at a concentration of **1M 200 pmol/ul**
- 3 For each nucleofection, prepare 6ml StemFlex + Revitacell or Rock inhibitor and coat 1× 6-well plate with geltrex/synthemax/matrigel for **01:00:00** at **37 °C**
- 4 For each nucleofection, prepare complete P3 solution by adding **20 µL** of Supplement to **90 µL** of P3 solition in a 1.5ml Eppendorf

B. Pre-assembly of Cas9 RNP

- 5 For each nucleofection, mix **4 µL** of sgRNA (**16 µg** total) and **2 µL** of recombinant Cas9 protein (**20 µg** total), in 1.5ml Eppendorf tube
- 6 Incubate at RT for **00:30:00** to **00:45:00** . *Note: begin harvesting iPSC after 15mins*
- 7 Immediately prior to nucleofection, add **1 µL** (**200 pmol**) of Alt-R HDR modified ssODN to the pre-assembled Cas9 RNP.

C. Delivery of Cas9 RNP and ssODN using Nucleofector

- 8 Switch on the nucleofector unit and select "X-unit" and single cuvette. Use the "Primary Cell P3" program and set the pulse code to **"CA-137"**
- 9 Collect the iPSC from an 80% confluent well with pre-warmed accutase. Count cells and pellet 8×10^5 cells by centrifuging at **300 x g, 20°C, 00:05:00** and aspirate supernatant.

- 10 While the cells are spinning, add the  1 μL of ssODN to the pre-assembled Cas9 RNP Eppendorf.
- 11 After pelleting cells and supernatant removed, gently resuspend the cell pellet in  100 μL complete P3 Solution and transfer the cells to the Eppendorf containing Cas9 RNP+ssODN. Transfer the cell suspension to the nucleocuvette. Tap the cuvette on a flat surface to eliminate bubbles and nucleofect immediately using the "Primary Cell P3" program and "CA-137" pulse code.
- 12 Immediately after the nucleofection, transfer the cells into one well of pre-coated, warm 6-well plate containing 2-3 ml of Stemflex + revitacell/Rock. Inh.
- 13 Add HDR enhancer to a final concentration of  30 micromolar (μM) (1:100 from the 3mM stock solution). Culture the cells in a  32 $^{\circ}\text{C}$ incubator (cold shock for 2 days: important for homozygous gene editing).
- 14 Change media the next day to StemFlex without revitacell/rock inh and HDR enhancer
- 15 Return cells to  37 $^{\circ}\text{C}$ incubator on day 3

D. Plating single cells to isolated clones

- 16 Once the cells in the 6-well plate are 80% confluent, dissociate them with accutase into a single cell suspension, centrifuge at  210 x g, 20 $^{\circ}\text{C}$, 00:05:00 . Remove supernatant and resuspend cell pellet in StemFlex + revitacel/rock.inh
- 17 Count the cells and plate between 800 – 1000 single cells in geltrex/synthemax/matrigel-coated 10cm petri dishes containing 10ml Stemflex + revitacel/rock inh.
- 18 Transfer cells to  37 $^{\circ}\text{C}$ incubator.
- 19 Change the media the next day and daily with Stemflex to remove revitacell/rock inhibitor

E. Picking single cell-derived clones

- 20 With the aid of a dissecting microscope and a p200, individual colonies are picked and transferred into 2× 96-well plates, one for colonies expansion, the other for DNA lysis and genotyping.
- 21 For each 96 colonies to be picked, prepare 1x geltrex coated 96-well and 15ml of Stemflex+Rock inh/revitacell and warm to RT in the hood.
- 22 Use the p200 tip to break apart the colony from the plate and collect the cell fragments in a volume of  50 μL of media. Transfer each colony to a well of a U-bottom 96-well plate.
- 23 Once 96 colonies are picked, add  150 μL of Stemflex+rock inh/revitacell to each well of the U-bottom plate with a multichannel.
- 24 With the use of the multichannel set to 100ul, pipette cells up and down 10 times in the U-bottom plate. Transfer  100 μL of the cell suspension to geltrex coated 96-well plate (for colony expansion) and  100 μL to 96-well plate for lysis and DNA extraction.
- 25 The following day, replace the media in the expansion 96-well plate with fresh Stemflex without revitacell/rock inh.

F. Genomic DNA extraction from 96-well plates

- 26 For each 96-well plate to be lysed, add  10 mg of proteinase K powder (Sigma-Aldrich, P6556) to  10 mL of lysis buffer ( 50 millimolar (mM) KCl,  10 millimolar (mM) Tris-HCl pH 8,  2 millimolar (mM) MgCl_2 ,  0.45 % volume (v/v) NP40,  0.45 % volume (v/v) Tween20).
- 27 With a multichannel pipettor, remove media from confluent wells of the Synthemax-coated 96-well plate and wash each well once with  150 μL room temperature D-PBS.



- 28 Add  100 μL of lysis buffer containing 1 mg/ml proteinase K, seal the 96-well plate and incubate at  60 °C for at least  04:00:00 .
- 29 While the plate is still warm, transfer lysates to a 96-well PCR plate, seal the plate and heat inactivate the proteinase K by incubating the plate at  95 °C for  00:10:00 in a thermocycler.
- 30 While the samples are warm, prepare a 1:20 dilution of the lysates in water for setting up PCR reactions. Store the undiluted lysates at  -80 °C . The diluted samples can be kept at  4 °C for a few days.