



Mar 15, 2024

CRISPR gRNAs cloning

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v9yj3pg3e/v1>

Anita Adami¹

¹Laboratory of Molecular Neurogenetics, Department of Experimental Medical Science, Wallenberg Neuroscience Center and Lund Stem Cell Center, BMC A11, Lund University, 221 84 Lund, Sweden.



Danielle Daft

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v9yj3pg3e/v1>

Protocol Citation: Anita Adami 2024. CRISPR gRNAs cloning . protocols.io

<https://dx.doi.org/10.17504/protocols.io.x54v9yj3pg3e/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 23, 2022

Last Modified: May 31, 2024

Protocol Integer ID: 57294

Keywords: CRISPR gRNAs cloning, gRNA oligonucleotides, ASAPCRN, procedure of crispr grnas cloning, crispr grnas cloning, crispr, procedure, crispr grna

Funders Acknowledgements:

Aligning Science Across Parkinson's through the Michael J. Fox Foundation for Parkinson's Research

Grant ID: ASAP-000520

Swedish Research Council

Grant ID: 2018-02694

Swedish Brain Foundation

Grant ID: FO2019-0098

Cancerfonden

Grant ID: 190326

Barncancerfonden

Grant ID: PR2017-0053

NIHR Cambridge Biomedical Research Centre

Grant ID: NIHR203312

Swedish Society for Medical Research

Grant ID: S19-0100

National Institutes of Health

Grant ID: HG002385

Swedish Research Council

Grant ID: 2021-03494

Swedish Research Council

Grant ID: 2020-01660

Abstract

This protocol details the procedure of CRISPR gRNAs cloning.

Attachments



[gfccbg5jf.pdf](#)

40KB



gRNA oligonucleotides design

- 1 To design your gRNAs, use CRISPick portal

(<https://portals.broadinstitute.org/gppx/crispick/public>).

- 1.1 Order the Oligos with specific overhangs for BsmBI cloning.

- 1.2 Insert the designed 20bp target gRNA sequence between the overhangs.

Forward oligo: 5' CACCG.....20 bp target.....-3'

Reverse oligo: 5' AAAC.....20 bp.....C 3'

gRNA oligonucleotides cloning

- 2 **Preparation of the gRNA oligonucleotides.**

- 2.1 Spin oligonucleotide tubes briefly.



- 2.2 Dilute to 100 micromolar (μM) solution with water.

- 2.3 Vortex, leave for some minutes, and vortex again.



- 2.4 Annealing of oligonucleotides:



- 100 micromolar (μM) of oligo A (forward)
- 100 micromolar (μM) of oligo B (reverse)
- 2 μL 10x NEB buffer 2
- water up to 20 μL total reaction volume



2.5 Denature at 95 °C for 00:05:00 , then cool down slowly.

5m

Note

Recommended: Turn the heating block off and leave the tubes in it for 02:00:00 - 03:00:00

2.6 When they have reached Room temperature , spin down.



3 **Prepare the assembly reaction for each oligonucleotide in individual PCR tubes containing:**



1. 100 ng backbone of lentiviral plasmid of choice (make sure it includes the AmpR gene for selection)
2. 1 µL of annealed gRNA oligonucleotide from step 1
3. 1 µL BsmBI/Esp3I restriction enzyme. FAST DIGEST
4. 1 µL T4 DNA ligase
5. 2 µL 10x T4 ligase buffer (to a final concentration of 1x)
6. Nuclease-free water up to 20 µL total reaction volume

3.1 Incubate the reaction in a thermal cycler with the following conditions:

1h

- 10 cycles 00:05:00 at 37 °C .
 00:10:00 at 22 °C .
- Hold for 00:30:00 at 37 °C .
- Hold for 00:15:00 at 75 °C .
- Keep at 4 °C .



Plasmid transformation & preparation

4h 32m 45s



- 4 Thaw Stbl3 or homemade top10 competent bacteria On ice .
- 4.1 Add 2 μL of the ligation reaction from step 2 to the bacteria On ice .
- 4.2 Mix a little by tapping the tube carefully a couple of times.
- 4.3 Keep On ice for 00:30:00 . 30m
- 4.4 To transform, dip the tubes in a 42 °C water bath for exactly 00:00:45 . 45s
- 4.5 Put the tubes back On ice for 00:02:00 . 2m
- 4.6 Transfer the bacteria to 250 μL pre-warmed or Room temperature soc media in a ventilated 15 ml falcon tube.
- 4.7 Incubate at 37 °C with shake for 01:00:00 . 1h
- 4.8 Spread everything on pre-warmed ampicillin+ agar plates.
- 4.9 Incubate at 37 °C Overnight . 1h
- Note

Only successfully transformed colonies will grow on the plate.
- 5 The day after, pick up 3 different colonies for each of the plasmids from the agar plates and prepare 3 minipreps. Incubate the minipreps Overnight on shake at 37 °C . 1h



6 Isolate the plasmid from the minipreps using the GeneJet Plasmid Miniprep kit (ThermoFisher).

6.1 Measure the DNA concentration.

6.2 Digest the DNA using the AflIII (BspTI) restriction enzyme to linearize the plasmid.



Note

This restriction site is part of the LTR found in lentiviral plasmids.

6.3 Check for the correct plasmid size on 1% agarose gel.

6.4 If the plasmids have the correct size, send for sequencing 1 or 2 for each cloned gRNA..

7 If the sequencing confirms the correct plasmid sequence, use one of the sequenced miniprep for each gRNA to prepare a maxiprep. Incubate the maxipreps at 37 °C .

Overnight

1h



8 Isolate the plasmid from the maxipreps using the NucleoBond Xtra Midi Plus Ef (ThermoFisher).

8.1 Measure the DNA concentration.

8.2 Digest the DNA using the AflIII (BspTI) restriction enzyme to linearize the plasmid.



8.3 Check for correct plasmid size on 1% agarose gel, and send for sequencing.