

Nov 30, 2023

CRISPRa tiling screens

DOI

<https://dx.doi.org/10.17504/protocols.io.q26g7p9bqgwz/v1>

Sean R. McCutcheon¹, Adam M. Swartz¹, Michael C. Brown¹, Alejandro Barrera¹, Christian McRoberts Amador¹, Keith Siklenka¹, Lucas Humayun¹, Maria A. ter Weele¹, James M. Isaacs¹, Andrea Daniel¹, Timothy E. Reddy¹, Andrew Allen¹, Smita K. Nair¹, Scott J. Antonia¹, Charles A. Gersbach¹

¹Duke University

Andrea Daniel: This protocol was adapted from the work of Sean McCutcheon and colleagues in the Gersbach lab at Duke University.

Gersbach Lab



Andrea R Daniel

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g7p9bqgwz/v1>

Protocol Citation: Sean R. McCutcheon, Adam M. Swartz, Michael C. Brown, Alejandro Barrera, Christian McRoberts Amador, Keith Siklenka, Lucas Humayun, Maria A. ter Weele, James M. Isaacs, Andrea Daniel, Timothy E. Reddy, Andrew Allen, Smita K. Nair, Scott J. Antonia, Charles A. Gersbach 2023. CRISPRa tiling screens. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.q26g7p9bqgwz/v1>

Manuscript citation:

McCutcheon, S.R., Swartz, A.M., Brown, M.C. *et al.* Transcriptional and epigenetic regulators of human CD8⁺ T cell function identified through orthogonal CRISPR screens. *Nat Genet* (2023). <https://doi.org/10.1038/s41588-023-01554-0>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 30, 2023

Last Modified: November 30, 2023

Protocol Integer ID: 91655

Keywords: CRISPRa, dSaCas9, Jurkat cells, IL2RA, tiling screen, tiling guide rna screen, guide rna screens in jurkat cell, activity of dsacas9, rna, dsacas9, using promoter, jurkat cell, cell

Funders Acknowledgements:

NIH

Grant ID: HG012053

Abstract

This protocols describes methods for characterizing the activity of dSaCas9 as a activator using promoter tiling guide RNA screens in Jurkat cells.

Construction of CRISPRa Jurkat lines

- 1 Polyclonal dSaCas9^{VP64} and ^{VP64}dSaCas9^{VP64} Jurkat cell lines were generated by transducing Jurkat cells with lentivirus encoding for either dSaCas9^{VP64}-2A-PuroR or ^{VP64}dSaCas9^{VP64}-2A-PuroR.
- 2 Cells were selected for 5 days using 0.5 µg ml⁻¹ of puromycin.

IL2RA CRISPRa tiling screen

- 3 After selection, 1×10^6 dSaCas9^{VP64} and ^{VP64}dSaCas9^{VP64} Jurkat cells were plated and transduced in triplicate with the *IL2RA* gRNA library lentivirus at a low multiplicity of infection (MOI).
- 4 Cells were expanded for 10 days, selected for Thy1.1 using a CD90.1 Positive Selection Kit (StemCell Technologies), and stained for Thy1.1 and IL2RA.
- 5 Transduced cells in the lower and upper 10% tails of IL2RA expression were sorted for subsequent gRNA library construction and sequencing.
- 6 All replicates were maintained and sorted at a minimum of 500× coverage.

gRNA sequencing

- 7 Genomic DNA was isolated using Qiagen's DNeasy Blood and Tissue Kit. Genomic DNA was split across 100 µl PCR reactions (25 cycles at 98 °C for 10 s, 60 °C for 30 s, and 72 °C for 20 s) with Q5 2× Master Mix and up to 1 µg of genomic DNA per reaction.
- 8 PCRs were pooled together for each sample and purified using double-sided (SPRI) bead selection at 0.6× and 1.8×.
- 9 Libraries were run on a High Sensitivity D1000 tape (Agilent) to confirm amplicon size and quantified using Qubit's dsDNA High Sensitivity assay.
- 10 Libraries were diluted to 2 nM, pooled together at equal volumes, and sequenced using Illumina's MiSeq Reagent Kit v2 (50 cycles).



- 11 Primers are available in Supplementary Table 5 of McCutcheon et al. Nature Genetics, 2023. <https://doi.org/10.1038/s41588-023-01554-0>

Processing gRNA sequencing and gRNA analysis

- 12 FASTQ files were aligned to custom indexes for each gRNA library (generated from the bowtie2-build function) using Bowtie 2 (ref. [67](#)).
- 13 Counts for each gRNA were extracted and used for further analysis in R.
- 14 Individual gRNA enrichment was determined using the DESeq2 (ref. [68](#)) package to compare gRNA abundance between groups for each screen.

Protocol references

67. Lagmead, B. & Salzberg, S. L. Fast gapped-read alignment with Bowtie 2. Nat. Methods 9, 357–359 (2012).

68. Love, M. I., Huber, W. & Anders, S. Moderated estimation of fold change and dispersion for RNA-seq data with DESeq2. Genome Biol. 15, 550 (2014).