

Jan 14, 2020

Cryopreservation

DOI

<https://dx.doi.org/10.17504/protocols.io.8rihv4e>

Michael Eadon¹, Daria Barwinska¹, Ying-Hua Cheng¹, Michael J. Ferkowicz¹, Samir V. Parikh², Brad H. Rovin², John P. Shapiro², Pierre C. Dagher¹, Tarek M. El-Achkar¹

¹Indiana University School of Medicine; ²Ohio State University Wexner Medical Center

KPMP

Tech. support email: info@kpmp.org



Michael Eadon

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8rihv4e>

Protocol Citation: Michael Eadon, Daria Barwinska, Ying-Hua Cheng, Michael J. Ferkowicz, Samir V. Parikh, Brad H. Rovin, John P. Shapiro, Pierre C. Dagher, Tarek M. El-Achkar 2020. Cryopreservation. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.8rihv4e>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 26, 2019

Last Modified: April 11, 2020

Protocol Integer ID: 29194

Keywords: cryopreservation samples of kidney tissue, kpmp biopsy tissue, kpmp pilot nephrectomy, cryopreservation sample, kidney tissue, other tissue interrogation site, other tissue interrogation sites for further interrogation, μm sections for imaging study, nephrectomy, iu biopsy, biopsy, shared tissue block, imaging data, kidney, tissue block, proteomic, imaging, imaging study, tissue, efficient use of the edge tissue

Abstract

Samples of kidney tissues are cryopreserved, sectioned and processed for various analyses. KPMP biopsy tissue undergoes initial processing (e.g. OCT embedding) at the Recruitment Sites, prior to distribution to Tissue Interrogation Sites. KPMP pilot nephrectomies were cryopreserved at University of Michigan. IU biopsies and nephrectomies were cryopreserved as described below. The tissue is sectioned according to a predetermined order: Bulk, LMD, miRNA, proteomics at 12 μm and then 50 μm sections for imaging studies as described in [Section 4.4](#). This allows efficient use of the edge tissue and maximization of imaging data that will anchor the LMD information with the remainder of the biopsy to be sent to other Tissue Interrogation Sites for further interrogation. Shared tissue blocks received from other Tissue Interrogation Sites will be sectioned in the reverse order.

Guidelines

This protocol is to be used for 0.5-10 mm^3 pieces of kidney tissue, including local nephrectomies and biopsies.

Cryopreservation

1

Note

This protocol is to be used for 0.5-10mm³ pieces of kidney tissue, including local nephrectomies and biopsies.

Kidney tissues are acclimated to OCT for  00:03:00 and then transferred to a cryomold with partially frozen OCT in the bottom (on a block of dry ice).

2 The tissue is completely covered in OCT, avoiding bubbles.

3 When the OCT is completely frozen, the tissue block is wrapped in parafilm (aiming to expel all air), then placed in a sealed bag and stored at  -80 °C

Cryosectioning Procedure

4

Note

This protocol is to be used with kidney tissue preserved in OCT and store at  -80 °C

All work is performed in a manner that limits RNA contamination and necessitates use of clean disposable gloves and face mask, as well as ensuring the cleanliness of all surfaces (RNase Away, Ambion, Cat #10328011).

5 1.2 µm Leica PPS-membrane slides (Leica, Cat# 11505268) used for LMD are exposed to UV light (in a tissue culture laminar flow hood) for 30 minutes, immediately prior to cryosectioning. Slides are stored at room-temp for optimal tissue adherence.

6 Cryostat is cooled to  -22 °C . The work surfaces are cleaned and a new cutting blade is installed.

- 7 A small slide box (cleaned with RNase Away) is placed inside the cryostat chamber to store slides with freshly cut tissue.
- 8 Specimen in OCT is adhered to a tissue holder and allowed to equilibrate for a few minutes to reach the chamber temperature and strengthen the adhesion between the OCT block and the holder. This process is aided via use of a heat extractor.
- 9 The specimen is cut at the following thickness:
 - a) 12 μm (2 section) – placed in Eppendorf tube for Bulk analysis
 - b) 12 μm (8 sections) – affixed to the specialized Leica LMD slide; one nephrectomy section per slide or two biopsy sections per slide. The slides are stored in $-80\text{ }^{\circ}\text{C}$, with a desiccant cartridge (Bel-Art, Cat# F42046-0000). To prevent any moisture from accumulating inside the slide box, it is further stored in a tightly closed Ziplock® bag.
 - c) 12 μm (1 section) – affixed to a glass slide for Periodic acid-Shiff staining
 - d) 12 μm (2 sections) – miRNA
 - e) 12 μm (6 sections) – For proteomics, tissue is affixed to specialized PEN-membrane LMD slides; one nephrectomy section per slide or two biopsy sections per slide. The slides are stored in $-80\text{ }^{\circ}\text{C}$, with a desiccant cartridge prior to shipment to OSU.
 - f) 50 μm (2 sections) – stored in tissue culture plate with 4% PFA for 24hr, followed by transfer of tissue into a scintillation vial with 0.25% PFA for long term storage. The tissue in 0.25% PFA is stored at 4°Celsius at all times when not in use. All sections per donor can be placed into the same storage container. Subsequent imaging of 50 μm slices is described in section 4.4
- 10 The LMD membrane slide adapter is used to assist with tissue adherence and collection.
- 11 Each slide is labeled with a specimen ID, date, and slide number.
- 12 Slides are used within 10 days from the date of cryosectioning.
- 13 For specimens transferred elsewhere for processing, the tissue is shipped (Mon-Thurs) overnight on dry ice to be received at the destination facility the following business day.