


Jun 01, 2023

Cryptococcus neoformans DNA Extraction Method

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l2522jl1y/v1>

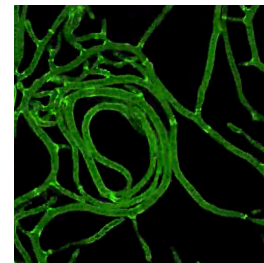

Amy Gladfelter¹

¹Duke University



Amy Gladfelter

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l2522jl1y/v1>

Protocol Citation: Amy Gladfelter 2023. Cryptococcus neoformans DNA Extraction Method. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.4r3l2522jl1y/v1>

Manuscript citation:

Based on Pitkin et al. (1996) Microbiology 142: 1557-1565. Results in ~0.5-2 mg DNA.

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 02, 2018

Last Modified: June 01, 2023

Protocol Integer ID: 17443

Keywords: cryptococcus neoformans dna extraction method extraction method, genomic dna from cryptococcus neoforman, cryptococcus neoforman, genomic dna, extraction, dna

Abstract

Extraction method to obtain genomic DNA from Cryptococcus neoformans.

Attachments



[DNAextraction_Southe...](#)

51KB

Guidelines

Extraction buffer (100 ml)

	Stock Solution	Add	Final Concentration
	1 M Tris-HCl, pH 7.5	10 ml	100 mM
	5 M NaCl	14 ml	0.7 M
	0.5 M EDTA	2ml	10 mM
	CTAB powder	1g	1%
	B-mercaptoethanol (14 M)	1ml	1%
	Water	73 ml	

CTAB is mixed alkyltrimethyl ammonium bromide, **Sigma cat.# M7635**. This takes time to go into solution. You can also use solid NaCl rather than a 5 M solution if that is easier. The buffer lasts four- six months at room temperature. The buffer seems to work better for long strands and spooling if β -mercaptoethanol is added just prior to use.

Safety warnings

 See SDS (Safety Data Sheet) for hazards and safety guidelines.



1 Grow a  50 mL YPD culture overnight ( 16:00:00), shaking at  30 °C .

2 Pellet cells in tabletop centrifuge in a 50 ml disposable tube.

Note

Optional: wash pellet with water and repeat spin.

3 Freeze cells at  -20 °C to  -80 °C for <30 min, then dry in a freeze drying machine.

4 Add the equivalent of  3 mL to  5 mL of 2 mm glass beads and vortex/shake until the cell pellet is broken and a fine powder is created.

5 **In fume hood**, add  10 mL CTAB extraction buffer (see Guidelines) and mix.

6 Incubate at  65 °C for  00:30:00 .

7 **In fume hood**, add  10 mL chloroform and gently mix for approximately  00:01:00 .

8 Spin in a table top centrifuge for  00:10:00 (2,500 – 3,000 rpm).

9 Remove supernatant (c. 7 ml) and add to an equal volume of isopropanol in a 15 ml disposable tube.

10 Gently rock back and forward to mix.



If the DNA precipitates in strands and clumps, spool out with a glass pipette and transfer to eppendorf containing  1 mL 70% ethanol.

Otherwise, spin in a table top centrifuge for  00:10:00 , pour off supernatant and use  1 mL 70% ethanol to wash DNA pellet and transfer it to an eppendorf tube.

11 Spin sample in microcentrifuge for 5-10 minutes. Remove ethanol and allow to air dry.

12 Resuspend DNA in either water or TE buffer (c. 500 μ l).

RNase can be added to final concentration of 20 μ g/ml if needed.

Run  1 μ L on an agarose gel to check concentration and quality.