

Jan 27, 2026

CTAB–Chloroform DNA Extraction from Ethanol-Preserved Freshwater Phytobentos

DOI

<https://dx.doi.org/10.17504/protocols.io.dm6gpzxwdlzp/v1>

Pavla Urbankova¹, Jan Štěpka¹

¹Faculty of Science, J. E. Purkyně University in Ústí nad Labem, Pasteurova 3632/15, 40096 Ústí nad Labem, Czech Republic



Pavla Urbankova

Univerzita Jana Evangelisty Purkyně

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.dm6gpzxwdlzp/v1>

Protocol Citation: Pavla Urbankova, Jan Štěpka 2026. CTAB–Chloroform DNA Extraction from Ethanol-Preserved Freshwater Phytobentos. protocols.io <https://dx.doi.org/10.17504/protocols.io.dm6gpzxwdlzp/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: July 20, 2024

Last Modified: January 27, 2026

Protocol Integer ID: 103761

Keywords: metabarcoding, CTAB, DNA extraction, chloroform, isopropanol, chloroform dna extraction from ethanol, extraction of edna, chloroform dna extraction, freshwater phytobentos this protocol, single chloroform extraction, preserved freshwater phytobento, stream water sample, based extraction, preserved river, ethanol, ctab

Abstract

This protocol describes a CTAB-based extraction of eDNA from particulate material pelleted from an ethanol-preserved river/stream water sample. It uses a single chloroform extraction, isopropanol precipitation, two 70% ethanol washes, and elution in Qiagen EB buffer. The protocol is adapted from Strand et al. (2019) with modifications for a pellet-based workflow.

Guidelines

- Contamination prevention (recommended): Use filter tips, wear gloves, change gloves frequently, and decontaminate work surfaces with a bleach-based solution before and during the work.
- Pellet handling: DNA pellets may be small and can be nearly invisible to the naked eye. Use extra care when removing the isopropanol supernatant after precipitation and during the 70% ethanol wash steps to avoid pellet loss.
- Optional clarification step: If the recovered aqueous phase remains visibly turbid after chloroform separation, an additional chloroform extraction can be performed to improve phase clarity before precipitation.



Materials

Reagents

- CTAB buffer (pH 8.0): 100 mM
- Tris-HCl, 20 mM EDTA, 1.4 M NaCl, 2% (w/v) CTAB (no additives)
- Proteinase K (20 mg/mL)
- Chloroform, pure (p.a.)
- Isopropanol (p.a.)
- 70% ethanol (prepared from p.a. ethanol + autoclaved deionized water)
- Qiagen EB buffer (10 mM Tris-Cl, pH ~8.5)

Consumables

- 2.0 mL microcentrifuge tubes (standard snap-cap "Eppendorf" type)
- Filter tips
- Gloves

Equipment

- Microcentrifuge for 2.0 mL tubes (able to run at ~13,500 rpm / max speed)
- Temperature-controlled device (heat block, water bath, or incubator) capable of 65 °C
- Vortex mixer
- Vacuum concentrator (optional; alternative is heat block drying)

Safety warnings

- !
 - Chloroform is toxic and volatile: handle in a fume hood; dispose of organic waste properly.
 - Ethanol and isopropanol are flammable: keep away from ignition sources.
 - Bleach-based disinfectants are irritants/corrosive: wear gloves and avoid mixing with acids.

Before start

- Pre-cool isopropanol and 70% ethanol at -20 °C.
- Preheat a heat block, water bath, or a temperature-controlled incubator (suitable for microcentrifuge tubes) to 65 °C.
- Decontaminate the work surface and equipment and organise a clean workflow to minimise contamination.



Pellet preparation from ethanol-preserved sample (2 mL aliquot)

33m 30s

- 1 Resuspend particulates in the ethanol-preserved sample container by manual shaking until homogenized. 30s
- 2 Transfer  2 mL of homogenized ethanol mixture into a 2.0 mL microcentrifuge tube. 30s
- 3 Centrifuge at  13500 rpm, Room temperature, 00:20:00 (max speed). 21m
- 4 Carefully decant the supernatant without disturbing the pellet. 30s
- 5 Dry the pellet for  00:10:00 using either a vacuum concentrator or a heat block (open cap), ensuring ethanol is removed. 11m

CTAB lysis and Proteinase K digestion

2h 35m

- 6 Add  500 μ L CTAB buffer (room temperature) to the dried pellet. 30s
- 7 Vortex continuously for  00:15:00 . 16m
- 8 Freeze at  -20 $^{\circ}$ C until fully frozen (typically  01:00:00); samples may be stored at  -20 $^{\circ}$ C  Overnight and processed the next day. 1h 1m
- 9 Thaw at  65 $^{\circ}$ C for  00:15:00 . 16m
- 10 Add  4 μ L 20 μ L Proteinase K ( 20 mg/mL) and vortex briefly. 30s
- 11 Incubate at  65 $^{\circ}$ C for  01:00:00 . 1h 1m



Chloroform extraction

19m

- 12 Add  400 μL chloroform and mix gently with the pipette tip. 1m
- 13 Centrifuge at  13500 rpm, Room temperature, 00:15:00 . 16m
- 14 Carefully transfer ~  300 μL of the upper aqueous phase to a new 2.0 or 1,5 mL tube using a fresh filter tip, avoiding the interphase. 2m

Isopropanol DNA Precipitation

33m

- 15 Add  400 μL ice-cold isopropanol (stored at  -20 °C) and mix by inverting the tube several times.
- 16 Incubate  00:15:00 at  4 °C . 16m
- 17 Centrifuge for  13500 rpm, Room temperature, 00:15:00 . 16m
- 18 Remove supernatant carefully without losing the pellet. 1m

Ethanol washes

15m

- 19 Add  500 μL ice-cold 70% ethanol (stored at  -20 °C) and briefly vortex. 1m
- 20 Centrifuge at  13500 rpm, Room temperature, 00:05:00 . 6m
- 21 Carefully remove supernatant. 30s



22 Repeat ethanol wash once more  [go to step #19](#) .

7m 30s

Drying and elution

12m

23 Dry the pellet (open cap) for ~  00:10:00 in a vacuum centrifuge or on a  65 °C heat block (avoid airborne contamination).

11m

24 Dissolve DNA pellet in  30 µL EB buffer (or sterile milliQ water), vortex and spin down. Let the DNA dissolve for at least 1 hour before further analysis (or store in fridge/freeze).

1m

Protocol references

Strand, D. A., Johnsen, S. I., Rusch, J. C., Agersnap, S., Larsen, W. B., Knudsen, S. W., Møller, P. R., & Vrålstad, T. (2019). *Monitoring a Norwegian freshwater crayfish tragedy: eDNA snapshots of invasion, infection and extinction*. *Journal of Applied Ecology*, 56(7), 1661–1673. <https://doi.org/10.1111/1365-2664.13404>