



Jun 09, 2018

Version 2

CTAB genomic DNA extraction from Arabidopsis leaf material V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.quidwue>

Pip Wilson¹, Diep Ganguly², Xin Hou¹, Barry Pogson¹

¹Australian National University; ²University of Pennsylvania

Pogson Group



Diep R R Ganguly

The Australian National University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.quidwue>

Protocol Citation: Pip Wilson, Diep Ganguly, Xin Hou, Barry Pogson 2018. CTAB genomic DNA extraction from Arabidopsis leaf material. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.quidwue>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 09, 2018

Last Modified: June 09, 2018

Protocol Integer ID: 12906

Keywords: genomic DNA, Arabidopsis, dna extraction from arabidopsis leaf material extraction, arabidopsis leaf material extraction, extraction of genomic dna, dna extraction, arabidopsis leaf material, genomic dna, dna, ctab, genomic, extraction

Abstract

Extraction of genomic DNA from Arabidopsis leaf material.

Guidelines

Gives reasonable quality and yield of gDNA, typically would use for PCRs, sequencing, and cloning but not for next-generation sequencing.

Materials

MATERIALS

✕ RNase A **Qiagen Catalog #19101**

✕ EDTA

✕ 1.5 mL Eppendorf tubes

✕ water

✕ Ethanol

✕ NaCl **Merck MilliporeSigma (Sigma-Aldrich) Catalog #53014**

✕ Hexadecyltrimethylammonium bromide **Merck MilliporeSigma (Sigma-Aldrich) Catalog #H6269**

✕ Tris-HCl (Tris-Hydrochloride), 100gm **Promega Catalog #H5121**

✕ 2-Propanol (IsoPropanol) **Bio Basic Inc. Catalog #PC8601.SIZE.4L**

✕ Tris-EDTA, pH 8.0 **Ambion Catalog #AM9849**

✕ Chloroform **Merck MilliporeSigma (Sigma-Aldrich) Catalog #366919-1L**

✕ Centrifuge

✕ Water bath set to 65°C

Safety warnings

! Perform chloroform steps in fume hood.



Before start

Make CTAB buffer:

- 2% (w/v) CTAB
- 1.4 M NaCl
- 0.1 M Tris-HLC pH 8

Grind leaf tissue into fine powder using mortar and pestle or Qiagen tissue lyser (place 1/8" steel ball bearing into tube with tissue sample).



Pre-heat 2% CTAB buffer at 65 °C in water bath

- 1 Heat only desired volume for use (~300 µl/100mg leaf tissue)

Add 300 µl CTAB buffer to each sample and mix well (vortex)

- 2
 - Adjust CTAB volume = 300 µl/100mg tissue
 - Can add 3 µl RNase A solution (100 µg/µl) to each sample if desired.

Incubate samples at 65°C (water bath) for at least 30 mins with occasional mixing by inversion of tubes.

- 3 Can be up to several hours or as needed based on input tissue and desired yields.

Remove samples from water bath and allow to cool to room temp.

- 4
 - Can place on ice/in fridge to speed up.

Add 300 µl chloroform to each sample. Mix well (vortex or shaking).

- 5
 - Perform in fume hood.

Centrifuge 5 - 15 min at 20,000g and transfer upper aqueous layer to new tube.

- 6
 - Transfer ~100 - 200 µl (lower volume = less chance of contamination from organic layer).
 - Depending on gDNA use, yield/cleanliness required, can repeat steps 5-6.

Add equal volume ice cold 2-propanol and incubate for 30 min @ -20 °C

- 7
 - Can incubate longer or at -80 °C depending on yield required.



Centrifuge @ 20,000g for 5 - 15 min to form a pellet and discard supernatant using pipette.

8

Add 500 μ l 70 - 80 % Ethanol to sample and centrifuge @ 9,000 - 10,000g for 5 min. Pipette off ethanol (as much as possible) and air dry for 1-2 min.

9

Resuspend gDNA in desired volume of water or Tris-EDTA buffer. Determine concentration by Nanodrop or Qubit or run on agarose gel to check quality.

10