



May 24, 2023

Culture and transfection of iPSC-derived neurons for live-imaging of axonal cargoes

 [Cell reports](#)

DOI

dx.doi.org/10.17504/protocols.io.x54v9dj4zg3e/v1

Dan Dou^{1,2}, Alexander Boecker³, Erika L.F. Holzbaur^{1,2}

¹Department of Physiology, Perelman School of Medicine, University of Pennsylvania, Philadelphia, PA 19104, USA;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, USA;

³Department of Neurology, University Medical Center Goettingen, 37077 Goettingen, Germany



Dan Dou

University of Pennsylvania

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

[Create free account](#)

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v9dj4zg3e/v1>

External link: <https://doi.org/10.1016/j.celrep.2023.112448>

Protocol Citation: Dan Dou, Alexander Boecker, Erika L.F. Holzbaur 2023. Culture and transfection of iPSC-derived neurons for live-imaging of axonal cargoes. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.x54v9dj4zg3e/v1>

Manuscript citation:

Boecker, C.A., and Holzbaur, E.L.F. (2021). Hyperactive LRRK2 kinase impairs the trafficking of axonal autophagosomes. *Autophagy* 00, 1–3. Boecker, C.A., Olenick, M.A., Gallagher, E.R., Ward, M.E., and Holzbaur, E.L.F. (2020). ToolBox: Live Imaging of intracellular organelle transport in induced pluripotent stem cell-derived neurons. *Traffic* 21, 138–155. Fernandopulle, M.S., Prestil, R., Grunseich, C., Wang, C., Gan, L., and Ward, M.E. (2018). Transcription Factor-Mediated Differentiation of Human iPSCs into Neurons. *Curr. Protoc. Cell Biol.* 79, e51.

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 13, 2022

Last Modified: June 01, 2024

Protocol Integer ID: 71294

Keywords: iPSC, iNeuron, live-imaging, axon, ASAPCRN, imaging of axonal cargo, glutamatergic neuron, excitatory glutamatergic neuron, excitatory glutamatergic neurons for the purpose, transport of axonal cargo, neuron differentiation from human ipsc, derived neuron, axonal cargo, differentiation of neuron, transfection of ipsc, disk confocal microscopy, neuron, transfect human ipsc, confocal microscopy, derived excitatory, microscopy, ipscs for differentiation, human ipsc, imaging, ipsc, live imaging

Funders Acknowledgements:

ASAP

Grant ID: ASAP-000350

Abstract

Here, we plate, culture, and transfect human iPSC-derived excitatory glutamatergic neurons for the purpose of observing transport of axonal cargoes under spinning disk confocal microscopy. Protocol is largely as previously described (Boecker et al., 2020, 2021; Fernandopulle et al., 2018). For preceding differentiation of neurons, see "Protocol: Piggybac-mediated stable expression of NGN2 in iPSCs for differentiation into excitatory glutamatergic neurons" and "Protocol: iNeuron differentiation from human iPSCs."

Attachments



[550-1146.pdf](#)

168KB

Materials

Materials

Equipment	
35 mm Dish No. 1.5 Coverslip 20 mm Glass Diameter Uncoated	NAME
Coverslip	TYPE
Mattek	BRAND
p35g-1-5-20-c	SKU
https://www.mattek.com/store/p35g-1-5-20-c-case/	LINK

Reagents

- PLO (CATALOG)
- Borate buffer (CATALOG)
- BrainPhys media (CATALOG)
- NT-3 (CATALOG)
- BDNF (CATALOG)
- B-27 supplement (CATALOG)
- Mouse laminin (CATALOG)
- 5-Fluoro-2'-deoxyuridine
- Uridine

Troubleshooting



Culture and transfection of iPSCderived neurons for live-imaging of axonal cargoes

30m

- 1 In advance, prepare 10x PLO stock.



	A	B
	PLO	50 mg
	0.1M borate buffer	50 mL

Note

Store 10x PLO stock at -80 °C .

- 2 The day before plating, coat imaging dishes with 1x PLO solution (10x PLO stock diluted in ddH₂O).



Note

It is only necessary to fully coat the glass center of the imaging dish.

- 3 The day of plating, remove PLO solution from imaging dishes and wash twice with ddH₂O.



- 3.1 Add 2 mL of iNeuron culture media.



BrainPhys supplemented with

	A	B
	BDNF	10 ng/mL
	NT-3	10 ng/mL
	Laminin	1 µg/mL




	A	B
	B-27 supplement	1x

- 3.2 Place dishes in cell culture incubator for >  00:30:00 .

30m



- 4 Rapidly thaw cryopreserved iNeurons in  37 °C water bath.

Note

Retrieve vial to tissue culture hood when only a small amount of ice remains visible.

- 5 Centrifuge to remove freezing media and resuspend cell pellet in iNeuron culture media.



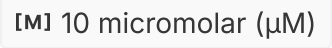
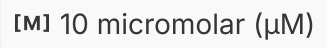
BrainPhys supplemented with

	A	B
	BDNF	10 ng/mL
	NT-3	10 ng/mL
	Laminin	1 µg/mL
	B-27 supplement	1x

- 6 Count cells and plate 300k neurons per 35 mm imaging dish.



- 6.1 Add cells dropwise to the center area of the dish (so that they sink onto the glass, PLO-coated center).


- 7 For Piggybac-delivered NGN2 neurons, include  10 micromolar (µM) 5-Fluoro-2'-deoxyuridine and  10 micromolar (µM) uridine at the time of plating to prevent survival of mitotic cells.

**Note**

These drugs were removed 24 hours after plating.



- 8 Store neurons in cell culture incubator. Perform partial change of iNeuron media twice per week.



- 9 On DIV18 (~  72:00:00 prior to imaging), transfect iNeurons for imaging.

3d

**Note**

Transfection conditions may require optimization, but a typical transfection will use  4 μL Lipofectamine Stem and  1 μg of plasmid DNA. Plasmids with the PGK or EF1 α promoters express best in iNeurons.