

Jun 09, 2025

CycloneSEQ long-read RNA-seq Library Preparation and Sequencing Protocol

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.e6nvwq3o7vmk/v1>

Qunfei Guo¹, youliang Pan¹, Wei Dai¹, Wei Jiang¹

¹BGI

GigaScience Press

BGI



shi shuaizhen

BGI research WhHan



Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvwq3o7vmk/v1>

Protocol Citation: Qunfei Guo, youliang Pan, Wei Dai, Wei Jiang 2025. CycloneSEQ long-read RNA-seq Library Preparation and Sequencing Protocol. protocols.io <https://dx.doi.org/10.17504/protocols.io.e6nvwq3o7vmk/v1>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 07, 2025

Last Modified: June 09, 2025

Protocol Integer ID: 219762

Keywords: CycloneSeq, Long-read Sequencing, RNA-seq, Genomics, Transcriptomics, animal rna on the cycloneseq platform, cycloneseq platform, cycloneseq, length transcriptome library preparation, read rna, extracted animal rna, transcriptome, rna, seq library preparation, sequencing protocol, sequencing, sequencing protocol this protocol, seq

Abstract

This protocol is designed for full-length transcriptome library preparation and sequencing of extracted animal RNA on the CycloneSEQ platform.



Genomic DNA Removal

- 1 Treat 100–500 ng total RNA with 5× gDNA Digester Mix in a nuclease-free tube.
 - Incubate at 42°C for 2 min.
 - Centrifuge briefly to collect liquid.

First-Strand cDNA Synthesis

- 2 Prepare 20 µL reaction mix containing:
 - 4× Hifair SuperMix
 - Random Primers (50 µM)
 - Treated RNA template

Incubate at 72°C for 5 min.

- Heat-inactivate at 85°C for 5 sec.

Note: For high-GC RNA templates, increase incubation temperature to 60°C.

CycloneSEQ Library Preparation

- 3 Process cDNA into sequencing libraries using the CycloneSEQ Universal Library Prep Set:
 - DNA Repair (end-repair/dA-tailing)
 - Adapter Ligation (CycloneSEQ-specific adapters)

Follow the standard whole-genome sequencing workflow steps provided in the kit.

Library Quality Control

- 4 Quantify libraries using a Qubit fluorometer.

Assess fragment size distribution (e.g., using Agilent TapeStation).

Sequencing

- 5 Load libraries onto the CycloneSEQ WuTong02 platform.

Run sequencing using manufacturer-recommended parameters (e.g., 48-hr run, >Q20 accuracy).