



Feb 13, 2024

Cylinder test in rats

DOI

<https://dx.doi.org/10.17504/protocols.io.kxygx36zdg8j/v1>

eduard.bentea¹, María Sanchiz Calvo¹, Veerle Baekelandt¹

¹Katholieke Universiteit Leuven



Eduard Bentea

KU Leuven

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kxygx36zdg8j/v1>

Protocol Citation: eduard.bentea , María Sanchiz Calvo, Veerle Baekelandt 2024. Cylinder test in rats. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.kxygx36zdg8j/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 13, 2024



Last Modified: June 01, 2024

Protocol Integer ID: 95156

Keywords: ASAPCRN, cylinder test in rat, cylinder test in rats protocol, cylinder test, nigrostriatal dopaminergic pathway, test, rats protocol, animal models with unilateral lesion, rat, asymmetry in forelimb use

Funders Acknowledgements:

ASAP (Aligning Science Across Parkinson's)

Abstract

Protocol for performing the cylinder test in rats. The cylinder test evaluates asymmetry in forelimb use, and can be used in animal models with unilateral lesions of the nigrostriatal dopaminergic pathway.



Test

1h

- 1 Bring cages to the behavioral room for at least  01:00:00 prior to the test for habituation  1h
- 2 Place each rat in a glass cylinder (20-cm wide), surrounded by mirrors to allow a full 360 degree view. Videotape from the front for  00:05:00 or until the rat performs a minimum of 20 weight-bearing forepaw contacts with the cylinder walls.  5m
- 3 Perform blinded analysis from the recorded videos of the number of contralateral forepaw contacts, and express these as a percentage of total forepaw contacts.