

Jun 03, 2020 Version 5

DAB Detection of Biocytin Labeled Tissue V.5

DOI

dx.doi.org/10.17504/protocols.io.bg5y jy7w

Allen Institute for Brain Science¹

¹Allen Institute

BICCN / BICAN

Allen Institute for Brain S...



Allen Institute

Allen Institute

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.bg5y jy7w

Protocol Citation: Allen Institute for Brain Science 2020. DAB Detection of Biocytin Labeled Tissue. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bg5y jy7w>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 03, 2020

Last Modified: March 17, 2021

Protocol Integer ID: 37784

Keywords: DAB, biocytin, diaminobenzidine, staining, coverslipping, mounting, PF0285,



Abstract

This protocol describes the process for diaminobenzidine (DAB) detection of biocytin filled cells. This protocol is optimized for use with brain slices cut at 350 μm thick, in which cells are first filled with biocytin (i.e., post-electrophysiological recording), fixed in 4% PFA/2.5% glutaraldehyde, and transferred to PBS until ready to stain.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



PF0285_DAB_Detection.



311KB

