



May 15, 2024

Version 6

DAB Detection of Biocytin Labeled Tissue V.6

DOI

dx.doi.org/10.17504/protocols.io.eq2ly3xkegx9/v6

Allen Institute for Brain Science¹

¹Allen Institute

Allen Institute for Brain Science

Tech. support

[Click here to message tech. support](#)



Allen Institute

Allen Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.eq2ly3xkegx9/v6>

Protocol Citation: Allen Institute for Brain Science 2024. DAB Detection of Biocytin Labeled Tissue. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.eq2ly3xkegx9/v6> Version created by **Allen Institute**

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 15, 2024

Last Modified: May 15, 2024

Protocol Integer ID: 99893

Keywords: DAB, biocytin, diaminobenzidine, staining, coverslipping, mounting, PF0285, dab detection of biocytin, detection of biocytin, dab detection, process for diaminobenzidine, use with brain slice, brain slice, tissue this protocol, dab, electrophysiological recording, filled cell, national institutes of health, cell,

Abstract

This protocol describes the process for diaminobenzidine (DAB) detection of biocytin filled cells. This protocol is optimized for use with brain slices cut at 350 μm thick, in which cells are first filled with biocytin (i.e., post-electrophysiological recording), fixed in 4% PFA/2.5% glutaraldehyde, and transferred to PBS until ready to stain.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



PF0285_DAB_Detection.



222KB

Troubleshooting

