


Mar 31, 2020

## DAB Staining of FFPE Slides

DOI

[dx.doi.org/10.17504/protocols.io.bdiei4be](https://dx.doi.org/10.17504/protocols.io.bdiei4be)

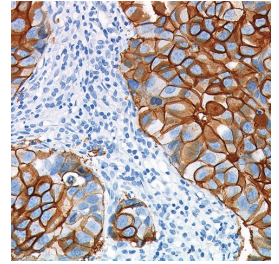

Jerelyn Nick<sup>1</sup>, Franchesca Farris<sup>1</sup>, Marda Jorgensen<sup>1</sup>

<sup>1</sup>University of Florida

Human BioMolecular Atlas Program (HuBMAP) Method Development Community  
Tech. support email: [Jeff.spraggins@vanderbilt.edu](mailto:Jeff.spraggins@vanderbilt.edu)



Marda Jorgensen

OPEN  ACCESS



DOI: [dx.doi.org/10.17504/protocols.io.bdiei4be](https://dx.doi.org/10.17504/protocols.io.bdiei4be)

**Protocol Citation:** Jerelyn Nick, Franchesca Farris, Marda Jorgensen 2020. DAB Staining of FFPE Slides. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.bdiei4be>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** March 11, 2020

**Last Modified:** October 02, 2020

**Protocol Integer ID:** 34086

**Keywords:** IHC, DAB, antibody validation, HubMAP, FFPE tissue ,



## Abstract

Standard IHC staining, such as this DAB protocol, is included in the validation of new antibodies at the University of Florida Tissue Mapping Center, TMC-Florida/Zurich. This procedure includes all of the reagents necessary to perform a DAB based immunohistochemistry reaction. DAB Substrate (3,3'-diaminobenzidine) produces a brown reaction product in the presence of peroxidase (HRP) enzyme.

## Attachments



New ELITE HRP Run

Re...

16KB

## Guidelines








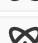











Microwave retrieval conditions are specific to the microwave in use. Correct temperature is achieved when small bubble form along the plastic coplin, but a boil does not occur.

A water jacket (beaker of room temperature water at a level that is 1cm from the cap level of the retrieval solution plastic coplin jar) is used to equilibrate the temperature and avoid hot spots during the retrieval.

Do not allow slides to become dry during staining procedure.

## Materials

### MATERIALS

-  Water
-  Ethanol, 200 proof
-  Hydrogen Peroxide, 30% **Fisher Scientific Catalog #H325-500**
-  Methanol **Fisher Scientific Catalog #A452-4**
-  Xylene **Fisher Scientific Catalog #X3P-1GAL**
-  Coplin Staining Jar, Plastic Coplin staining jar, Holds ten slides back-to-back **Thermo Fisher Catalog #107**
-  Shandon®; Shandon-Mount®; **Thermo Fisher Catalog #9990435**
-  Hematoxylin OPTIK DK **Catalog #RS4576-A**
-  Bluing Reagent OPTIK **Catalog #RS-4363-B**
-  Aqueous Clarifier **Catalog # RS4361-B**
-  Vector DAB Staining Kit **Vector Laboratories Catalog #SK-4100**
-  Vector Elite ABC Kit **Catalog #PK-4000**
-  Immunohistochemistry Slide Humidity Chamber **Newcomer Supply Catalog #68432A**
-  Avidin/Biotin Blocking Kit **Vector Laboratories Catalog #SP-2001**
-  Normal Serum for Blocking (see guide below) **Vector Laboratories**
-  Biotinylated anti-Primary Species Secondary (See selection guide below) **Vector Laboratories**
-  Tween 20 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #P9416**
-  TRIS-buffered saline (TBS 10X) pH 7.6 **Fisher Scientific Catalog #AAJ62662K3**
-  Citrate Buffer pH 6.0 10X **Abcam Catalog #ab64214**

### SECONDARY ANTIBODY AND BLOCKING SERUM SELECTION INFORMATION:

Primary Antibody Species		Biotinylated Secondary Antibody		Blocking Serum
RABBIT		Goat anti-Rabbit (BA-1000)		Goat (S-1000)
MOUSE		Horse anti-Mouse (BA-2000)		Horse (S-2000)
RAT		Rabbit anti-Rat (BA-4000)		Rabbit (S-5000)
GOAT		Rabbit anti-Goat (BA-5000)		Rabbit (S-5000)
CHICKEN		Goat anti-Chicken (BA-9010)		Goat (S-1000)

All Catalog Numbrers Provided are for Vector Laboratories Products



It is easier to work with Tween-20 if a 10% solution is prepared in advance of performing this protocol.

## Safety warnings

❗ Xylene is flammable and a neurotoxin. Use in a fumehood, keep containers tightly closed when not in use, and avoid open flames.

Hydrogen peroxide is ractive and can cause skin/eye damage. Handle with appropriate protective gear. Collect waste for chemical disposal

Methanol is flammable and an aquatic life toxin . Avoid flames, dispose of as chemical waste, collect first rinsate as waste.

Handle heated retrieval solutions with hot hands or heat resistant gloves.

DAB is a suspected carcinogen. Appropriate care should be exercised when using this reagent including gloves, eye protection, lab coats, and good laboratory procedures. Dispose in accordance with local regulations.

## Before start

Replenish chemicals in slide deparaffinization/rehydration portion of staining line.

Prepare humidity chamber for slide incubations.

A bright field bench microscope is needed to monitor DAB development reaction.

Blocking serum and biotinylated secondary antibody must be compatible with primary antibody source species.



## Day 1 Prepare Slides for staining

1d

1 De-paraffinize and rehydrate slides.

Place slides in a xylene-compatible rack.

1.1 Immerse slide holder in first xylene station for 5 minutes. ⌚ 00:05:00

1.2 Immerse slide holder in second xylene station for 5 minutes. ⌚ 00:05:00

1.3 Immerse slide holder in first 100% Ethanol for 5 minutes. ⌚ 00:05:00

1.4 Immerse slide holder in second 100% Ethanol for 5 minutes. ⌚ 00:05:00

1.5 Transfer slide holder into 3% H<sub>2</sub>O<sub>2</sub> in methanol for 10 minutes. ⌚ 00:10:00

### Note

Prepare the 3% H<sub>2</sub>O<sub>2</sub> in methanol solution fresh each day (25ml of 30% H<sub>2</sub>O<sub>2</sub> added to 225ml of methanol).

1.6 Immerse slide holder in 95% Ethanol for 5 minutes. ⌚ 00:05:00

1.7 Immerse slide holder in 75% Ethanol for 5 minutes. ⌚ 00:05:00

1.8 Immerse slide holder in 50% Ethanol for 5 minutes. ⌚ 00:05:00

1.9 Immerse slide holder in deionized water for 5 minutes. ⌚ 00:05:00

1.10 Immerse slide holder in deionized water for second wash for 5 minutes. ⌚ 00:05:00



## Antigen Retrieval

1d

- 2 Perform heat mediated antigen retrieval using 1X citra buffer.

### Note

Alternative antigen retrieval methods can be used at this step based on individual antibody product sheet or laboratory experience/preferencve.

- 2.1 Remove slides from slide holder and place into a plastic Coplin Jar filled with 10mM pH6.0 Na citrate buffer.  
Slides may be placed back to back in Coplin jar.  
Citrate buffer level must be over the top edge of the slides.
- 2.2 Place the Coplin Jar in a 400ml beaker filled with water leaving a centimeter of space between water level and cap bottom edge. Leave the cap loose, but held in place so that steam will not remove it.
- 2.3 Place beaker / Coplin jar assembly into microwave oven and heat for 7 minutes at 50% power.
- 2.4 Remove the beaker and jar from the microwave and incubate slides in the citrate buffer on the bench top for 18 more minutes.  
During this incubation, make 1X TBST buffer by combing 90ml de-ionized water, 10ml 10X TBS, pH 7.6, and 1ml 10% Tween-20.
- 3 Remove slides one at a time from the Coplin Jar and dip in de-ionized water to remove residual citra buffer.
- 3.1 Place slides in a humidity chamber and immediately cover sections with 1X TBST buffer. Equillibrate for 5 minutes

## Blocking and Applying Primary

1d

- 4 Avidin block  
To make the avidin solution: combine 1mL of 1X TBST buffer with 6 drops of Avidin from vector avidin biotin kit.
- 4.1 Take the slide, tap off buffer and cover section with Avidin Solution for 20 minutes.


00:20:00



## 5 Biotin block

To make the biotin solution: combine 1mL of 1X TBST buffer plus 6 drops of Biotin from a vector avidin biotin kit.

### 5.1 Take the slide, tap off buffer and cover section with the Biotin solution for 20 minutes.

 00:20:00

### 5.2 Tap off biotin solution and cover sections in 1X TBST buffer .


## 6 Serum block

To make the serum block, combine 1mL of 1X TBST buffer and 20uL of Normal Serum .

### Note

The blocking solution is prepared with from the serum from the same animal species in which the biotinylated secondary antibody is made.

### 6.1 Tap off buffer and remove the excess moisture with gauze. Cover section with serum block and incubate for 60 minutes in the humidity chamber

 01:00:00

### 6.2 Remove excess serum with gauze.

## 7 Primary antibody incubation

### Note

Prepare desired dilutions of primary antibody in Antibody Diluent (Thermo Fisher)

### 7.1 Apply primary antibody solution directly over the section, cover with a plastic slip cut from a clean plastic bag surface and place in humidity chamber. Incubate in a 4°C



refrigerator overnight. Overnight

## Day 2 Detect Antibody Staining

1d

- 8 Remove chamber from the refrigerator, carefully lift off and dispose of plastic bag coverslips.  
Wash slides with 1X TBST buffer for 5 minutes at room temperature . While washing, prepare ABC Reagent and secondary antibodies.

### Note

ABC Reagent needs to be made at least 30 minutes before it is used. To make the solution, use 20uL of reagent A and 1mL of buffer. Mix, then add 20uL of reagent B. Mix again.

To make secondary solution combine 1ml of 1X TBST buffer with 15ul of serum matching species secondary is raised in. Add 5ul of Vector biotinylated secondary antibody. Secondary antibody must be directed against species of origin for primary antibody.

- 9 Incubation with biotinylated secondary antibody(s).

- 9.1 Tip buffer of slides and remove excess moisture with gauze. Cover sections with the secondary antibody solution and incubate for 30 minutes at room temperature in a humidity chamber. 00:30:00

- 10 Wash slide with 1X TBST buffer for 5 minutes. 00:05:00  
Tip buffer of of slides and remove excess moisture with gauze.

- 11 Cover the sections with ABC reagent (step 8) and incubate for 30 minutes at room temperature. 00:30:00

- 12 Wash slides in 1X TBST buffer for 5 minutes. 00:05:00  
Slides can be held in buffer until ready for DAB staining.

- 13 DAB reaction  
To make the DAB solution: add 1 drop of DAB to 1mL of Impact Diluent.

- 13.1 Wipe excess buffer from one slide. Place the slide on to the stage of a bright field microscope, and identify an area for observing DAB development.





- 13.2 Cover the section on the slide with DAB solution and incubate until desired intensity develops. Once desired intensity is reached, rinse and leave in water until ready to counterstain.

#### Note

Collect DAB waste for hazardous chemical disposal.

## Counterstain and mount

1d

- 14 Counterstain.  
Prepare staining line in keeping with the steps below. Remove slides from water and place them into a xylene compatible slide holder.
- 14.1 Dip slide holder into hematoxylin for 10 seconds. ⌚ 00:00:10
- 14.2 Rinse with running tap water by immersing slide holder into water until it is clear.
- 14.3 Immerse slide holder into Clarifier for 20 seconds. ⌚ 00:00:20
- 14.4 Immerse slide holder into water for 1 minute.
- 14.5 Immerse slide holder into Bluing Reagent for 1 minute. ⌚ 00:01:00
- 14.6 Immerse slide holder into water for 1 minute.
- 14.7 Immerse slide holder into 50% Ethanol for 5 minutes. ⌚ 00:05:00
- 14.8 Immerse slide holder into 75% Ethanol for 5 minutes. ⌚ 00:05:00



- 14.9 Immerse slide holder into 95% Ethanol for 5 minutes. ⌚ 00:05:00
- 14.10 Immerse slide holder into 100% Ethanol for 5 minutes. Transfer slide holder into a second 100% Ethanol staining station for another 5 minutes. ⌚ 00:10:00
- 14.11 Immerse slide holder into first xylene for 3 minutes. Transfer slide holder into second xylene station for another 3 minutes. ⌚ 00:06:00
- 15 Mount a coverslip on to the slide using a permanent mountant.

#### Note

Working in the fume hood, remove one side at a time from the Xylene and mount coverslip with approximately 2 drops of Shandon Mount. Blot the edge of the coverslipped slide on a paper towel to remove excess mountant and allow to dry in the fume hood.

- 16 When slides are dry, image using a brightfield microscopy system. The antibody staining will appear brown and nuclei will be counterstained light blue.