

Oct 07, 2020

Data Processing of Nano-DESI IMS data

DOI

<https://dx.doi.org/10.17504/protocols.io.bm2sk8ee>

yang1832¹, Julia Laskin¹

¹Purdue University

NanoDESI MSI Julia Las...



yang1832

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bm2sk8ee>

Protocol Citation: yang1832 , Julia Laskin 2020. Data Processing of Nano-DESI IMS data. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bm2sk8ee>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: October 05, 2020

Last Modified: October 07, 2020

Protocol Integer ID: 42802

Keywords: mass spectrometry imaging data, desi ims data scope, mass spectrometry, final imzml file, data processing, nano, data, imaging, im

Abstract

Scope:

Process a mass spectrometry imaging data into a final imzML file.



- 1 Import the raw data folder into PeakByPeak Imaging software under "Process Raw folder"
- 2 Set "Number of Processes" to 10.
- 3 Choose "File Format" as imzML.
- 4 Run.