

Dec 01, 2023

Version 2

Deepwell Reusual Protocol V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlko2x6v5r/v2>

Morris Jack¹

¹UoL



Morris Jack

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlko2x6v5r/v2>

Protocol Citation: Morris Jack 2023. Deepwell Reusual Protocol . [protocols.io](https://dx.doi.org/10.17504/protocols.io.j8nlko2x6v5r/v2)

<https://dx.doi.org/10.17504/protocols.io.j8nlko2x6v5r/v2>Version created by [Morris Jack](#)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 01, 2023



Last Modified: December 01, 2023

Protocol Integer ID: 91694

Keywords: deepwell, reusual protocol, protocol

Abstract

This is a protocol to reuse deepwells for sustainability practices

Growing cultures in a deep well

- 1 Cover a deep well with two layers of aluminium foil and tape the front and back to hold it securely. The tape also turns the aluminium foil into a lid for the deep well by temporarily peeling off the front tape.
- 2 Autoclave at 120C for 20min. After autoclaving, keep in a heated oven (> 50C) overnight for drying and desiccation.
- 3 Do not fill the deep wells more than 60% of their max volume if they are to be placed in a shaking incubator to avoid spillage.

Deep well Decontamination

- 4 Tape around the deep well over the aluminium foil to prevent spillage.
- 5 While carrying the deep well outside the lab space, enclose it in a waterproof autoclave bag. Keep sufficient absorbent paper on hand in case of accidental spillage.
- 6 In the autoclave facility, autoclave the deep wells at 120C for >30min to decontaminate.
- 7 Once autoclaved, the deep wells will be brought back to the lab
- 8 Carefully peel the aluminium foil and dispose the cultures and aluminium foil into a disposal bag containing paper roll to absorb the liquid.
- 9 Send decontaminated culture for incineration (yellow bin route).
- 10 Rinse the deep well plate with water and a brush to clean before future use.
- 11 Keep in a heated oven (> 50C) overnight for drying and desiccation