

Oct 22, 2025

Denaturing On-bead Digestion for LC-MS/MS

 [Nature Communications](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l6zp91gqe/v1>

Katja Piltti¹, Clinton Yu¹

¹University of California, Irvine



Katja Piltti

University of California, Irvine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l6zp91gqe/v1>

Protocol Citation: Katja Piltti, Clinton Yu 2025. Denaturing On-bead Digestion for LC-MS/MS. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bp2l6zp91gqe/v1>

Manuscript citation:

Piltti KM et al. C1q drives neural stem cell quiescence by regulating cell cycle and metabolism through BAI1. Nature Communications (In press)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 08, 2025

Last Modified: October 22, 2025

Protocol Integer ID: 229372

Keywords: On-bead digestion, Denaturing protein digestion, Pull-down assay, Protein purification, Magnetic bead pull-down, Affinity purification, NHS-SS-Biotin labeling, DSP crosslinker, Streptavidin magnetic beads, Biotinylation-based protein capture, Crosslinked protein complexes, Proteomics sample prep, LC-MS/MS preparation, bead digestion for protein purification, bead digestion for lc, protein purification, protein complexes for unbiased forward screen, bead digestion, using magnetic bead, interacting intracellular protein, compatible with the cell membrane, magnetic bead, using nanolc, protein, biotin, protein complex, cell membrane

Funders Acknowledgements:

Wings for Life

Grant ID: WFL-US-01/18

Abstract

A denaturing on-bead digestion for protein purification after pull-down using magnetic beads. The protocol is compatible with the cell membrane-permeable Ez-link NHS-SS-Biotin (ThermoScientific, PI21441) labeling and cell membrane-permeable DSP (ThermoScientific, PI22586). Step-by-step instructions for the pull-down are available in "Co-Immunoprecipitation/Pull-Down of C1q-Interacting Intracellular Proteins/Protein Complexes for Unbiased Forward Screen Using NanoLC-MS/MS"-protocol.

Guidelines

1. Wash beads with HPLC grade H₂O and remove the supernatant.
2. Add 1 bed volume of 2M urea with 25 mM ammonium bicarbonate in water to the magnetic beads with protein.
Note: Bed volume refers to the volume of the resin or magnetic beads. For example, if you have 100 μ L of a 50% slurry, the bed volume is 50 μ L once settled.
3. Add Dithiothreitol (DTT) to a final concentration of 2 mM and incubate for 15 minutes at 37°C.
4. Add iodoacetamide to a final concentration of 4 mM and incubate for 30 minutes in the dark at room temperature.
5. Add DTT to a final concentration of 2 mM and incubate for 15 minutes at 37°C to quench.
6. Add 1 bed volume of 25 mM ammonium bicarbonate to dilute the urea to a final concentration of 1M.
7. Add trypsin to a final concentration of 20 ng/ μ L. Incubate at 37°C for 8 hours to overnight. *Dissolved in 50 mM acetic acid.
8. Add formic acid to 0.1% to stop the digestion.
9. Collect the supernatant from the beads for a c18 cleanup.



Materials

- HPLC grade H₂O (Fisher Scientific, Cat# W5-4)
- 2M urea (Fisher Scientific, Cat# BP169-500)
- Ammonium bicarbonate (Sigma-Aldrich, Cat# 09830)
- Dithiothreitol (DTT) (Fisher Scientific, Cat# BP172-25)
- Iodoacetamide (Sigma-Aldrich, Cat# I6125)
- Trypsin (Fisher Scientific, Cat# PRV5111)
- 50 mM acetic acid (Millipore-Sigma, Cat# 5330010050)
- Formic acid (Fisher Scientific, Cat# 60-006-17)

Proceed with the following steps for protein denaturation, reduction, and alkylation:

- 1 Wash beads with HPLC-grade H₂O (Fisher Scientific, Cat# W5-4) and remove the supernatant.
- 2 Add 1 bed volume of 2M urea (Fisher Scientific, Cat# BP169-500) with 25 mM ammonium bicarbonate (Sigma-Aldrich, Cat# 09830) in water to the magnetic beads with protein. Note: Bed volume refers to the volume of the resin or magnetic beads. For example, if you have 100 μ L of a 50% slurry, the bed volume is 50 μ L once settled.
- 3 Add Dithiothreitol (DTT) (Fisher Scientific, Cat# BP172-25) to a final concentration of 2 mM and incubate for 15 minutes at 37°C.
- 4 Add iodoacetamide (Sigma-Aldrich, Cat# I6125) to a final concentration of 4 mM and incubate for 30 minutes in the dark at room temperature.
- 5 Add DTT to a final concentration of 2 mM and incubate for 15 minutes at 37°C to quench.
- 6 Add 1 bed volume of 25 mM ammonium bicarbonate to dilute the urea to a final concentration of 1M.
- 7 Add trypsin* (Fisher Scientific, Cat# PRV5111) to a final concentration of 20 ng/ μ L. Incubate at 37°C for 8 hours to overnight. *Dissolved in 50 mM acetic acid (Millipore-Sigma, Cat# 5330010050).
- 8 Add formic acid (Fisher Scientific, Cat# 60-006-17) to 0.1% to stop the digestion.
- 9 Collect the supernatant from the beads for a C18 cleanup. See Sample Desalting Using OMIX C18 Tips for LC-MC-protocol for details.