

Jan 07, 2022

Detection of Cryptosporidium in stool sample by PCR-RFLP

DOI

<https://dx.doi.org/10.17504/protocols.io.bxaxpifn>

botchiesenyo¹

¹University of Ghana

Ayi and Pawlowic Lab C...



mcpawlowic

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bxaxpifn>

Protocol Citation: botchiesenyo 2022. Detection of Cryptosporidium in stool sample by PCR-RFLP. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bxaxpifn>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: August 11, 2021

Last Modified: January 07, 2022

Protocol Integer ID: 52279

Keywords: detection of cryptosporidium, identification of cryptosporidium spp, cryptosporidium spp, cryptosporidium, stool sample by pcr, environmental microbiology, stool sample, restriction fragment length polymorphism assay, microbiology, pcr, natural mineral water, rflp

Abstract

Nested PCR-RFLP adapted from Nichols, R.A.B., Campbell, B.M. and Smith, H.V., 2003. Identification of *Cryptosporidium spp.* oocysts in United Kingdom noncarbonated natural mineral waters and drinking waters by using a modified nested PCR-restriction fragment length polymorphism assay. *Applied and environmental microbiology*, 69(7), p.4183.

Materials

- GoTaq MasterMix from Promega
- Primers



Nested PCR

12h 45m

- 1 Extract whole genomic DNA from stool samples.

3h

Positive control DNA for *Cryptosporidium parvum* may be obtained from ATCC. Alternatively you can extract genomic DNA from wild type *Cryptosporidium parvum* purchased from Bunchgrass Farms (Deary, Idaho, USA), Waterborne Inc (New Orleans, Louisiana, USA), or Sterling Labs (Tucson, Arizona, USA).

- 2 First round of PCR to amplify 18SrRNA

4h

- 2.1 Combine the following in a PCR tube:

30m

5 µl of DNA (from positive control, experimental sample, or water for negative control)

0.1 µM Forward Primer ("out NDIAGF2": CAATTGGAGGGCAAGTCTGGTGCCAGC)

0.1 µM Reverse Primer ("out NDIAGR2": CCTTCCTATGTCTGGACCTGGTGAGT)

1x GoTaq Master Mix (Promega)

Ultrapure water to bring total volume to 25 microliters

- 2.2 Thermocycler program:

3h 30m

95 °C -5 minutes,

30 cycles:

94 °C -30 seconds

58 °C - 1 minute

72 °C - 30 seconds

- 3 Second round of PCR to amplify 18SrRNA

4h 30m

- 3.1 Combine the following in a PCR tube:

30m

1 µl of PCR product from Step 2

0.1 µM Forward Primer ("DIAGF": AAGCTCGTAGTTGGATTTCTG)

0.1 µM Reverse Primer ("DIAGR": TAAGGTGCTGAAGGAGTAAGG)

1x GoTaq Master Mix (Promega)

Ultrapure water to bring total volume to 25 microliters

- 3.2 Thermocycler program:

4h

95 °C -5 minutes,

45 cycles:

94 °C -30 seconds

58 °C - 1 minute

50 °C - 1 minute



- 4 Run 7 μ l of PCR product from Step 3 on a 2% agarose gel and visualise by staining with ethidium bromide.

45m

RFLP

4h 45m

- 5 Combine the following and incubate at 37 °C for 4 hours:

4h

5 μ l PCR product from step 4 above

20u Restriction enzyme (Ssp1, Vsp1, Ase1)

1x final concentration Restriction enzyme digestion buffer (corresponds to enzyme used)

1x final concentration BSA

- 6 Run digested PCR product from Step 5 on a 2% agarose gel and visualize by staining with ethidium bromide.

45m