


Sep 27, 2023

Version 2

Differentiation of RGC Induced Neurons (RGC-iNs) V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.14egn2pqzg5d/v2>

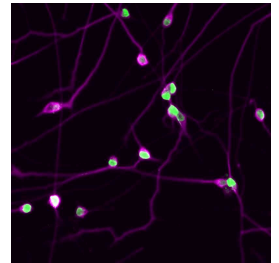

Devansh Agarwal^{1,2}, Kevin W. Mazo³, Karl Wahlin²

¹Department of Bioengineering, University of California San Diego, La Jolla, United States;

²Department of Ophthalmology, University of California San Diego, La Jolla, United States;

³Department of Biological Sciences, University of California San Diego, La Jolla, United States

Devansh Agarwal: Correspondence: d4agarwa@ucsd.edu;

Kevin W. Mazo: Correspondence: kewima99@gmail.com;

Karl Wahlin: Correspondence: kwahlin@ucsd.edu



Devansh Agarwal

Department of Bioengineering, University of California San D...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egn2pqzg5d/v2>

Protocol Citation: Devansh Agarwal, Kevin W. Mazo, Karl Wahlin 2023. Differentiation of RGC Induced Neurons (RGC-iNs).

protocols.io <https://dx.doi.org/10.17504/protocols.io.14egn2pqzg5d/v2> Version created by **Devansh Agarwal**

Manuscript citation:

Devansh Agarwal, Nicholas Dash, Kevin W. Mazo, Manan Chopra, Maria PA. Garcia, Amit Patel, Ryan M. Wong, Cairang Jia, Hope Do, Jie Cheng, Colette Chiang, Shawna L. Jurlina, Mike Perry, Jong Rho, Risa Broyer, Cassidy Lee, Robert N. Weinreb, Cezar Gavrilovici, Nick W. Oesch, Derek S. Welsbie, Karl J. Wahlin. Human retinal ganglion cell neurons generated by synchronous BMP inhibition and transcription factor mediated reprogramming. *npj Regen Med* **8**, 55 (2023). <https://doi.org/10.1038/s41536-023-00327-x>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 27, 2023

Last Modified: September 27, 2023

Protocol Integer ID: 88422

Keywords: differentiation of rgc induced neuron, rgc induced neuron, induced pluripotent stem cell, retinal ganglion cell, pluripotent stem cell, generating neuron, process of generating neuron, crispr, inducible polycistronic transcription factor gene cassette, containing human neurog2, neuron, human neurog2, cell, cas12a ribonucleoprotein

Funders Acknowledgements:

National Institutes of Health (NIH)

Grant ID: K99/R00EY024648

National Institutes of Health (NIH)

Grant ID: R01EY031318

National Institutes of Health (NIH)

Grant ID: R21EY031122

Abstract

This protocol is designed to convert human induced pluripotent stem cells (PSCs) into retinal ganglion cell induced neurons (RGC-iNs) using a doxycycline-inducible polycistronic transcription factor gene cassette containing human NEUROG2, ATOH7, ISL1, and POU4F2. The TetO-driven transgene cassette is integrated into the CLYBL safe harbor site using a CRISPR-Cas12a ribonucleoprotein. The process of generating neurons is greatly enhanced by the inclusion of the BMP blocker LDN-193189.

Guidelines

Apart from observation under the microscope, counting, and centrifugation, all steps should be carried out in a sterile biological safety cabinet.

Materials

Table: Key resources or reagents required.

A	B	C
REAGENT or RESOURCE	SOURCE	IDENTIFIER
40µm Cell Strainer, EASYstrainer (for generating single cell suspension during plating)	Greiner	Cat# 542040
Accutase (single cell passaging of hPSCs)	Sigma-Aldrich	Cat# A6964
B27 vitamin (neural supplement)	Thermo Fisher Scientific	Cat# 17504044
BDNF (growth factor for RGC growth and survival)	Qkine	Cat# Qk050
Blebbistatin (ROCK inhibitor for improving cell survival)	Sigma-Aldrich	Cat# B0560
BrainPhys Neuronal Medium (basal media for supporting long-term growth of neurons)	StemCell Technologies	Cat# 05790
CultureOne supplement (for enhancing neural conversion)	Thermo Fisher Scientific	Cat# A3320201
DMEM (basal media)	Thermo Fisher Scientific	Cat# 11965
DMEM/F12 50:50 (basal media)	Thermo Fisher Scientific	Cat# 11330
Doxycycline hyclate (antibiotic for transgene induction)	Sigma-Aldrich	Cat# D5207
F12 (basal media)	Thermo Fisher Scientific	Cat# 11765
GDNF (growth factor for enhancing RGC growth and neuronal survival)	Qkine	Cat# Qk051
Insulin-Human Recombinant (N2 supplement component)	Roche	Cat# 11376497001
L-ascorbic acid (N2 supplement component)	Sigma-Aldrich	Cat# A8960
LDN-193189 (pre-patterning BMP pathway inhibitor)	Sigma-Aldrich	Cat# SML0559
Matrigel® Growth Factor Reduced (GFR) Basement Membrane Matrix (cell attachment/differentiation of hPSCs)	Corning	Cat# 354230



	A	B	C
	mTeSR1 (maintenance and propagation of hPSCs)	Stem Cell Technologies	Cat# 85850
	N-2 Supplement (neural supplement)	Thermo Fisher Scientific	Cat# 17502048
	NEAA (non-essential amino acids for supporting neuronal growth)	Thermo Fisher Scientific	Cat# 11140
	Nicotinamide (NIC) (vitamin B3 supplement to enhance differentiation)	Sigma-Aldrich	Cat# 72340
	Poly-L-ornithine (PLO) hydrobromide (for neural attachment)	Sigma Aldrich	Cat# P3655
	Progesterone (N2 supplement component)	Sigma-Aldrich	Cat# P8783
	Putrescine dihydrochloride (N2 supplement component)	Sigma-Aldrich	Cat# P5780
	Sodium selenite (N2 supplement component)	Sigma-Aldrich	Cat# S5261
	Thiazovivin (alternate ROCK inhibitor for cell survival)	LC Labs	Cat# T-9753
	holo Transferrin Human (N2 supplement component)	Sigma-Aldrich	Cat# T0665
	GlutaMAX Supplement (auxiliary energy source for cells)	Thermo Fisher Scientific	Cat# 35050061

Before start

For media/reagent recipes see the last section of the protocol.

PSC expansion step

- 1 Grow PSCs in mTeSR1 under hypoxia ($[M] 5\%$ (v/v) O_2 / $[M] 10\%$ (v/v) CO_2) or normoxia ($[M] 20\%$ (v/v) O_2 / $[M] 5\%$ (v/v) CO_2) at $37\text{ }^{\circ}\text{C}$.

12-well plate (3.5 cm²): Plate 5,000 PSCs into each well of 12-well plates in mTeSR1 in the presence of $[M] 5\text{ }\mu\text{M}$ blebbistatin (blebb; 2,000x stock). Feed daily in mTeSR1 (without blebb) for ~4 more days. Typically, we get ~200,000 – 500,000 cells per well when cells are ready for passaging.

6-well plate (9.6 cm²): Plate 15,000 PSCs into each well of 6-well plates in mTeSR1 in the presence of blebb. Feed daily in mTeSR1 (without blebb) and grow for ~4 more days. Typically, we can get ~750,000 – 1,000,000 cells per well when cells are ready for passaging.

Note

For PSC expansion make sure that colonies are not overgrown (70-80 % confluent) and not touching.

Alternatively, $[M] 2\text{ }\mu\text{M}$ thiazovivin (10 mM or 5000x stock) can be used instead of blebb as a ROCK inhibitor.

Day -1: Priming of stem cells for neural induction

- 2 One day prior to neural induction, pre-coat TC plates overnight with $[M] 0.1\text{ mg/mL}$ poly-L-ornithine hydrobromide (PLO). To do this, add 1 mL $[M] 2\text{ mg/mL}$ PLO (20x) to 19 mL cell culture grade H₂O, use 1 mL per well for 6-well plates, 0.5 mL per well for 12 well plates and 0.25 mL per well for 24 well plates. Incubate overnight at $37\text{ }^{\circ}\text{C}$.

1d

Pre-treat stem cells by replacing mTeSR1 media with fresh mTeSR1 supplemented with $[M] 1\text{ }\mu\text{g/mL}$ doxycycline (dox; 1,000x stock) and $[M] 100\text{ nM}$ LDN-193189 (10,000x stock).

**Note**

For long-term experiments, you need better adhesion of cells so you can use a higher concentration of PLO. To do this, dilute  1 mL  2 mg/mL PLO (20x) into  9 mL cell culture grade H₂O for a final concentration of  0.2 mg/mL .

Day 0: Neural Induction

- 3 Wash overnight PLO-coated plates >3 times with culture grade H₂O. Let plates dry completely in the back of the TC hood for  01:00:00 , then coat with  1 % (v/v) Matrigel. For Matrigel coating, use  1 mL per well for 6-well plates,  0.5 mL per well for 12 well plates and  0.25 mL per well for 24 well plates. Incubate for >  03:00:00 at  37 °C .

4h

- 4 Prepare Neural Induction Medium (NIM) initiation cocktail with  2 µg/mL dox (500x stock),  100 nanomolar (nM) LDN (10,000x stock), 1x CultureOne (100x stock) and  5 micromolar (µM) blebb.

Note

Dox is light-sensitive, so keep aliquots cool after thawing and dark when not in use (up to ~1 month at  4 °C).

In addition, it is critical to prevent the drying of matrigel-coated dishes after the matrigel has been added.

- 5 Incubate the cells with prewarmed Accutase (~  00:05:00) in the incubator for  00:12:00 The volume of Accutase to use is 1/2 the volume that you maintain the cells in (e.g.,  1 mL per well of a 6 well plate,  0.5 mL per well of a 12 well plate and  0.25 mL per well of a 24 well plate).
- 6 Gently rinse the wells using a P1000 micropipette and pipet up and down 3 times to further break up the cell clumps into single cells.

17m



- 7 Put the cells into a conical  1.5 mL or  5 mL tube with 2 times the volume of prewarmed mTeSR1 with [M] 5 micromolar (μM) blebb (e.g.  1 mL Accutase +  2 mL mTeSR1) to quench the Accutase, then pellet the cells for  00:05:00 at  80 x g . 5m
- 8 Aspirate the supernatant and then resuspend the cell pellet in prewarmed NIM + [M] 5 micromolar (μM) blebb. 2m
- 9 Filter cells with a  40 μm cell strainer. Count the cells with a hemocytometer. 5m
- 10 Plate cells onto PLO/matrigel-coated plates. 5m
To make a 6-well plate (**9.6 cm²/well; 57.6 cm² total; 7,000 cells/cm²**): Add 403,200 cells into  12 mL (67,200 cells per well) of NIM initiation cocktail (NIM + dox, LDN, CultureOne, blebb) in a  15 mL conical tube, mix well and distribute across the wells ( 2 mL per well).
To make a 12-well plate (**3.5 cm²/well; 42 cm² total; 7,000 cells/cm²**): Add 294,000 cells into  12 mL (24,500 cells per well) of NIM initiation cocktail in a  15 mL conical tube, mix well and distribute across the wells ( 1 mL per well).
To make a 24-well plate (**1.9 cm²/well; 45.6 cm² total; 7,000 cells/cm²**): Add 319,200 cells into  12 mL (13,300 cells per well) of NIM initiation cocktail in a  15 mL conical tube, mix well and distribute across the wells ( 0.5 mL per well).

Day 1: Maintenance

- 11 Do nothing.

Day 2: Feed – add 1/3 of media

- 12 For 6 well plate: Add  1 mL **NIM +** [M] 1 $\mu\text{g}/\text{mL}$ **dox + 1xCultureOne** to each well. 5m
For 12 well plate: Add  0.5 mL **NIM + dox + CultureOne** to each well.
For 24 well plate: Add  0.25 mL **NIM + dox + CultureOne** to each well.



Note

Do this very carefully by adding media to the sides of the dish. If you are not very careful cells will detach.

Day 3: Maintenance

13 Do nothing.

Day 4: Feed – exchange 1/3 of media

- 14 For 6 well plate: Remove  1 mL media and replace with fresh  1 mL **NIM +**  1 µg/mL **dox + 1xCultureOne + NIC** ( 10 millimolar (mM)). 5m
- For 12 well plate: Remove  0.5 mL media and replace with fresh  0.5 mL **NIM + dox + CultureOne + NIC**.
- For 24 well plate: Remove  0.25 mL media and replace with fresh  0.25 mL **NIM + dox + CultureOne + NIC**.

Beyond day 4, plates need to be fed every other day.

Note

This is the last day to add CultureOne.

For all subsequent exchanges:

Neurons tend to easily dissociate from the dish, so be very careful when aspirating. Take care to aspirate and dissociate by tilting the dish so that the medium accumulates on one side. Then, aspirate/dispense with the pipette directed toward the wall of the dish (i.e., away from the cells at the bottom).

Day 6: Feed – exchange 1/3 of media

- 15 For 6 well plate: Remove  1 mL media and replace with fresh  1 mL **BrainPhys + B27 (1x) +**  1 µg/mL **dox + BDNF** ( 50 ng/mL) + **GDNF** ( 10 ng/mL) + **NIC** ( 10 millimolar (mM)). 5m

For 12 well plate: Remove  0.5 mL media and replace with fresh  0.5 mL

BrainPhys + B27 + dox + BDNF + GDNF + NIC.

For 24 well plate: Remove  0.25 mL media and replace with fresh  0.25 mL

BrainPhys + B27 + dox + BDNF + GDNF + NIC.

Note

This is the last day to add dox.

Day 8: Feed – exchange 1/3 of media

- 16 For 6 well plate: Remove  1 mL media and replace with fresh 1  1 mL **BrainPhys + B27 (1x) + BDNF ([IM] 50 ng/mL) + GDNF ([IM] 10 ng/mL) + NIC ([IM] 10 millimolar (mM)).**

5m

For 12 well plate: Remove  0.5 mL media and replace with fresh  0.5 mL

BrainPhys + B27 + BDNF + GDNF + NIC.

For 24 well plate: Remove  0.25 mL media and replace with fresh  0.25 mL

BrainPhys + B27 + BDNF + GDNF + NIC.

- 17 For long-term experiments >1 week, continue feeding by 1/3 media exchange every other day with **BrainPhys + B27 + BDNF + GDNF + NIC.**

5m

Media/Reagent Recipes

- 18 **Neural Induction medium (NIM medium):**

	A	B	C
	Component	Catalog #	Volume
	DMEM/F12, HEPES	Thermo Fisher Scientific # 11330	485 ml
	N2 supplement (100x)	Thermo Fisher Scientific # 17502048 or recipe below	5 ml
	NEAA (non-essential amino acids, 100x)	Thermo Fisher Scientific # 11140	5 ml



A	B	C
GlutaMAX supplement (100x)	Thermo Fisher Scientific # 35050061	5 ml
		Total 500 ml

19 **N2 supplement (100X) recipe** - If making in-house add the following:

A	B	C	D	E	F	G
Component	Catalog #	100 ml	150 ml	200 ml	250 ml	500 ml
Transferrin	Sigma-Aldrich # T0665	1 g	1.5 g	2 g	2.5 g	5.0 g
Insulin	Roche # 113764970 01	50 mg	75 mg	100 mg	125 mg	250 mg
Progesterone	Sigma-Aldrich # P8783	63 µg	94.5 µg	126 µg	157 µg	315 µg
Putrescine	Sigma-Aldrich # P5780	161 mg	241.5 mg	322 mg	402.5 mg	805 mg
Sodium Selenite	Sigma-Aldrich # S5261	50.2 µg	75.3 µg	100.4 µg	125.5 µg	251 µg
DMEM/F12	Thermo Fisher Scientific # 11330	to 100 ml	to 150 ml	to 200 ml	to 250 ml	to 500 ml

20 **Reagent Stock Dilutions:**

Recombinant Human BDNF Protein (Qkine, Cat# Qk050): 50 µg/mL stock in 10 millimolar (mM) HCl with 0.1 % (v/v) BSA; 1,000X; use at 50 ng/mL . Store aliquots in -80 °C .

Recombinant Human GDNF Protein (Qkine, Cat# Qk051): 10 µg/mL stock in cell culture grade H2O with 0.1 % (v/v) BSA; 1,000X; use at 10 ng/mL . Store



aliquots in -80 °C .

Doxycycline (dox) (Sigma-Aldrich, Cat# D5207): Make 1 mg/mL stock (working concentration is 0.5-2 µg/mL) in cell culture grade ddH₂O and filter sterilize; 1,000X; use at 1 µg/mL . Store 1 mL aliquots at -20 °C . Protect from light.

LDN-193189 hydrochloride (LDN) (Sigma Cat# SML0559-5MG): 1 millimolar (mM) (10,000x):

$g = \text{Molecular Weight (g/mol)} * \text{Molarity (M)} * \text{Volume (L)}$

$0.005 \text{ g} = (406.48 \text{ g/mol}) * (0.001 \text{ M}) * (\text{Volume})$

$0.005 \text{ g} = 0.41 \text{ g/L} * \text{Volume}$

Volume = 0.012 L or 5 mg LDN in 12 mL DMSO; Store aliquots in

-80 °C .

Matrigel (GF reduced) (1 % (v/v)) (Corning, Cat# 354230):

Thaw stock 10ml bottle overnight On ice (wet) before aliquoting. Always keep matrigel and tubes on ice and never let it come to room temperature or it will gel.

Make 200 µL aliquots. Store aliquots in -80 °C .

Add 200 µL matrigel to 24 mL ice-cold DMEM/F12 (~ 1 % (v/v) final).

Add 1 mL per well of 6-well plate, 0.5 mL per well of 12-well plate or

0.25 mL per well of 24-well plate.

Nicotinamide (NIC) (Sigma Cat# 72340): 1 Mass Percent (100x) stock solution (10 millimolar (mM) working solution). Soluble in water to ~ 1g/10ml.

$g = \text{Molecular Weight (g/mol)} * \text{Molarity (M)} * \text{Volume (L)}$; $g = (122.12\text{g/mol}) * (1\text{M}) * (0.05\text{L})$

= 6.11 g NIC in 50 mL of DMEM/F12 (or water); filter sterilize and store at

4 °C .

poly-L-ornithine hydrobromide (PLO) - (Sigma Cat# P3655-500MG; mol wt 30,000-70,000 Da): 2 mg/mL (20x) stock:

Add 500 mg PLO to 250 mL cell culture ddH₂O.

Make 1 mL aliquots. Store aliquots in -80 °C .

Blebbistatin (blebb) (10 millimolar (mM) or 2,000x stock, Sigma Cat# B0560-1MG):



$g = \text{Molecular Weight (g/mol)} * \text{Molarity (M)} * \text{Volume (L)}$

$0.001\text{ g} = (292.33\text{ g/mol}) * (0.01\text{ M}) * (\text{Volume})$

$0.001\text{ g} = 2.92\text{ g/L} * \text{Volume}$

$\text{Volume} = 0.000342\text{ L} = 0.342\text{ mL} = 342\text{ }\mu\text{L}$

Dissolve  1 mg blebbistatin into  342 μL DMSO, divide into  20 μL aliquots

(store at  $-80\text{ }^{\circ}\text{C}$). Working concentration is  5 micromolar (μM) .

Protocol references

1. Fernandopulle, M. S. *et al.* Transcription Factor-Mediated Differentiation of Human iPSCs into Neurons: Rapid differentiation of iPSCs into neurons. *Current Protocols in Cell Biology* **79**, e51 (2018).
2. Nehme, R. *et al.* Combining NGN2 Programming with Developmental Patterning Generates Human Excitatory Neurons with NMDAR-Mediated Synaptic Transmission. *Cell Reports* **23**, 2509–2523 (2018).