

Aug 18, 2020

Direct ELISA for investigating the binding of peroxidase-labeled anti-chicken IgY conjugate with avian immunoglobulins

DOI

<https://dx.doi.org/10.17504/protocols.io.bjxykppw>

Angel Justiz-Vaillant¹

¹University of the West Indies St. Augustine

University of the West In...

angel.vaillant@sta.uwi.e...



Angel A Justiz-Vaillant

University of the West Indies St. Augustine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bjxykppw>

Protocol Citation: Angel Justiz-Vaillant 2020. Direct ELISA for investigating the binding of peroxidase-labeled anti-chicken IgY conjugate with avian immunoglobulins. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.bjxykppw>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: August 18, 2020

Last Modified: August 18, 2020

Protocol Integer ID: 40664

Keywords: avian immunoglobulin, immunoglobulin, binding of peroxidase, peroxidase, igy conjugate, direct elisa, chicken, present in the egg, egg, avian species, igy

Abstract

The peroxidase-labeled anti-chicken IgY conjugate cross-reacts with many IgY present in the egg of many and diverse avian species.

References

1. Justiz Vaillant AA, Ramirez N, Cadiz A, Ferrer B, Akpaka P, et al. (2013) Separation and Reactivity of Avian Immunoglobulin Y. J Chromat Separation Techniq 4: 173. doi:10.4172/2157-7064.1000173

Materials

MATERIALS

⊗ Anti-Chicken IgY, HRP Conjugate, 300ul **Promega Catalog #G1351**

⊗ Nunc™ 96-Well Polystyrene Round Bottom Microwell Plates, V 96 well plate, Non-Treated, clear, without lid, Sterile **Thermo Fisher Catalog #260210**

- 1 This ELISA is used to study the interaction of anti-chicken IgY-HRP conjugate with diverse avian immunoglobulins.
- 2 The 96 well microtitre plate is coated overnight at 4°C with 1 µg/µl per well of purified avian immunoglobulins or 50 µl of water soluble fraction from egg yolks of avian species in carbonate-bicarbonate buffer pH 9.6.
- 3 Then plate is treated with bovine serum albumin solution and washed 4X with PBS-Tween.
- 4 Then 50 µl of peroxidase-labeled-anti-chicken IgY conjugate diluted 1:15000 in PBS-non-fat milk is added to each well and incubated for 1.30h at RT. After that the plate is washed 4X with PBS-Tween.
- 5 Pipette 50 µl of 3,3',5,5' - tetramethylbenzidine (TMB; Sigma-Aldrich) to each well.
- 6 The reaction is stopped with 50 µl of 3M H₂SO₄ solution.
- 7 The plate is visually assessed for the development of colour and read in a microplate reader at 450 nm.
- 8 A cut-off point should be calculated as the mean of the optical density of negative controls x 2 SD.