

Aug 18, 2020

Direct ELISA for investigating the binding of Protein-A to immunoglobulins.

DOI

<https://dx.doi.org/10.17504/protocols.io.bjxqkpmw>

Angel Justiz-Vaillant¹, Monica F. Smikle²

¹University of the West Indies St. Augustine; ²University of the West Indies. Mona Campus

University of the West In...

angel.vaillant@sta.uwi.e...



Angel A Justiz-Vaillant

University of the West Indies St. Augustine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bjxqkpmw>

Protocol Citation: Angel Justiz-Vaillant, Monica F. Smikle 2020. Direct ELISA for investigating the binding of Protein-A to immunoglobulins. . protocols.io <https://dx.doi.org/10.17504/protocols.io.bjxqkpmw>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 18, 2020

Last Modified: August 18, 2020

Protocol Integer ID: 40656

Keywords: immunoglobulin, binding of protein, direct elisa, protein, elisa

Materials

MATERIALS



Nunc™ 96-Well Polystyrene Round Bottom Microwell Plates, V 96 well plate, Non-Treated, clear, without lid, Sterile **Thermo Fisher Catalog #260210**



Staphylococcal Protein-A **Merck MilliporeSigma (Sigma-Aldrich)**

- 1 This ELISA is used to study the interaction of protein-A (SpA) with diverse immunoglobulins.
- 2 The 96 well microtitre plate is coated overnight at 4°C with 1 µg/µl per well of purified immunoglobulins or 50 µl of any animal sera in carbonate-bicarbonate buffer pH 9.6.
- 3 Then plate is treated with bovine serum albumin solution and washed 4X with PBS-Tween.
- 4 Then 50 µl of peroxidase-labeled-protein-A conjugate diluted 1:5000 in PBS-non-fat milk is added to each well and incubated for 1.30h at RT. After that the plate is washed 4X with PBS-Tween.
- 5 Pipette 50 µl of 3,3',5,5' - tetramethylbenzidine (TMB; Sigma-Aldrich) to each well.
- 6 The reaction is stopped with 50 µl of 3M H₂SO₄ solution.
- 7 The plate is visually assessed for the development of colour and read in a microplate reader at 450 nm.
- 8 A cut-off point should be calculated as the mean of the optical density of negative controls x 2 SD.