

Mar 27, 2023

Version 1

DNA - Ball Python DNA Amplification with TFEC primers V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl4jb78lo5/v1>



Jose Avila Cervantes¹

¹Lolita Genomics



Jose Avila Cervantes

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl4jb78lo5/v1>

Protocol Citation: Jose Avila Cervantes 2023. DNA - Ball Python DNA Amplification with TFEC primers. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.3byl4jb78lo5/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: October 18, 2022

Last Modified: March 27, 2023

Protocol Integer ID: 71489

Keywords: ball python dna amplification, tfec primers pcr amplification, piedball morph in ball python, piedball morph, pcr amplification, ball python, aactcagagcactccatgacc right primer, python regius, dna, genotype, primer, left primer, tfec, amplification

Abstract

PCR amplification with Tfec set Exon 5 to genotype Piedball morph in Ball python (Python regius)

LEFT PRIMER AACTCAGAGCACTCCATGACC

RIGHT PRIMER CAGGTGTGCCCCTTTCATAA

1 REAGENTS

- 10X PCR Buffer
- MgSO₄ [M] 20 millimolar (mM)
- Taq Polymerase 5U/ul
- Primer Tfec exon5 Left [M] 10 micromolar (μM)
- Primer Tfec exon5 Right [M] 10 micromolar (μM)
- DNTP's [M] 10 micromolar (μM)
- Ultra Pure Water
- 100bp ladder (100-2,000 bp)
- SYBR Safe DNA stain
- Loading dye
- Agarose

2 EQUIPMENT

- DNA LoBind tubes 1.5mL
- PCR 8-Strip tubes 0.2mL
- Micropipettes
- Thermal Cycler

3 MASTER MIX

30s



Below is the recommended volume for one sample. Adjust for the number of samples to process and add 10% to account for pipetting error. Thaw all reagents at room temperature before using them except the Taq DNA polymerase, which has to be kept on ice at all times. Premix all reagents and add the Taq DNA polymerase at the end, vortex thoroughly for  00:00:30 seconds. Aliquot the master mix in to each tube

 28 μL and then add  2 μL of DNA sample. The final volume for each reaction is

 30 μL .

	A	B
	Reagent	Volume (uL)
	10x PCR Buffer	3
	MgSO ₄	4.5
	Taq Polymerase	0.3

	A	B
	Primer F	1.5
	Primer R	1.5
	DNTP's	0.6
	Template-DNA	2
	Water	16.6

Table 1. Master Mix for PCR.

4 PCR CONDITIONS



	A	B	C	D
	STEP	TIME	TEMPERATURE (°C)	
	Initial Denaturation	3 min	94	
	Denaturation	15 sec	94	25 Cycles
	Annealing	30 sec	53.2	
	Extension	30 sec	74	
	Final Extension	2 min	74	

Table 2. Thermocycler program



5 AMPLICON QUALITY CHECK

35m



1. Prepare a 1% agarose gel with 1X SYBR Safe DNA stain.
2. Add  4 μL of 100bp ladder (100-2,000 bp) in the first well.
3. Premix  4 μL of sample and  1 μL of 10X loading dye or  3 μL of sample and  1 μL of 6X loading dye. Load this mixture in each well.
3. Run the samples for  00:35:00 at a 100V.
4. Visualize in a trans illuminator and take a photo of the gel.
5. The product should be a clear, single band around 200 bp.