

Mar 27, 2023

Version 1

DNA - Ball Python DNA Extraction from sheds V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzb344vx1/v1>



Jose Avila Cervantes¹

¹Lolita Genomics



Jose Avila Cervantes

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzb344vx1/v1>

Protocol Citation: Jose Avila Cervantes 2023. DNA - Ball Python DNA Extraction from sheds. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.rm7vzb344vx1/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: October 15, 2022

Last Modified: March 27, 2023

Protocol Integer ID: 71398

Keywords: ball python dna extraction, dna from ball python, ball python, dna, extraction, using phenol, python regius, isoamyl alcohol, dry shed

Abstract

Protocol to extract DNA from Ball Python (Python regius) dry sheds using Phenol:Chloroform:Isoamyl Alcohol



1 EQUIPMENT

- Dry Bath / Heated Block
- Microcentrifuge
- DNA LoBind tubes 1.5mL
- Micropipettes
- Assorted pipette tips

2 REAGENTS

- Lysis Buffer ([M] 10 millimolar (mM) Tris-base, [M] 100 millimolar (mM) EDTA, 2% SDS, [M] 5 Mass Percent , NaCl, $\text{pH } 8$)
- TE Buffer (EDTA [M] 1 millimolar (mM) , Tris-Cl [M] 10 millimolar (mM))
- Proteinase K ([M] 20 mg/mL)
- Phenol/Chloroform/Isoamyl Alcohol 25:24:1 (v/v)
- Ethanol 100 %
- Ethanol 70 % (Freshly prepared)
- Tris Acetate-EDTA (TAE) Buffer
- SYBR Safe DNA stain
- Agarose
- Loading Dye
- Wide Range Ladder (100-12,000 bp)

3 PROTEINASE K DIGESTION

30s

1. Cut a small piece of fresh tissue (3-5mm). Put in a 1,5mL microtube. Add  900 μL of lysis buffer. Add  9 μL Proteinase K ([M] 20 mg/mL) and vortex thoroughly for  00:00:30 seconds.



2. Incubate at  55 °C overnight .

4 PHENOL/CHLOROFORM/ISOAMYL EXTRACTION

5m 30s

1. Add  500 μL of Phenol/Chloroform/Isoamyl Alcohol 25:24:1 (v/v) and vortex thoroughly for  00:00:30 seconds.



2. Centrifuge at room temperature for  00:05:00 at  16000 x g .



3. Carefully remove the upper aqueous phase (~  500 μL) and transfer the layer to a fresh tube. Be sure not to carry over any Phenol during pipetting.

5 ETHANOL PRECIPITATION

1h 50m



1. Add  1000 μL of  100 % volume Ethanol, invert the tube and place the tube at  $-80\text{ }^{\circ}\text{C}$ for  01:00:00 or at  $-20\text{ }^{\circ}\text{C}$  Overnight .

2. Centrifuge the sample at  $4\text{ }^{\circ}\text{C}$ for  00:30:00 at  16000 x g to pellet the DNA.

3. Carefully remove the supernatant without disturbing the pellet.

4. Add  150 μL of  70 % volume ethanol.

5. Centrifuge the sample at  $4\text{ }^{\circ}\text{C}$ for  00:05:00 at  16000 x g .

6. Carefully remove the supernatant without disturbing the pellet.

7. Repeat  70 % volume ethanol wash once and remove as much of the remaining ethanol as possible.

8. Dry the DNA at  Room temperature for  00:05:00 to  00:10:00 .

9. Re-suspend the DNA pellet in  200 μL of TE Buffer.

6 DNA QUALITY CHECK

40m



1. Prepare a 1% agarose gel with 1X TAE buffer and 1X SYBR Safe DNA stain.

2. Add  4 μL Wide Range Ladder (100-12,000 bp) in the first well.

3. Premix  9 μL of sample and  1 μL of 10X loading dye or  5 μL of sample and  1 μL of 6X loading dye. Load this mixture in each well.



3. Run the samples for  00:40:00 at a 100V.
4. Visualize in a trans illuminator and take a photo of the gel.