

Nov 23, 2020

Version 1

## DNA extraction for human microbe samples. V.1

 [GigaByte](#)

 Peer-reviewed method

 In 8 collections

DOI

[dx.doi.org/10.17504/protocols.io.bcmriu56](https://dx.doi.org/10.17504/protocols.io.bcmriu56)

Lilan Hao<sup>1</sup>

<sup>1</sup>BGI-Shenzhen, Shenzhen 518083, China.

GigaScience Press

BGI



Lilan Hao

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bcmriu56>

External link: <https://doi.org/10.46471/gigabyte.9>



**Protocol Citation:** Lilan Hao 2020. DNA extraction for human microbe samples.. **protocols.io**

**<https://dx.doi.org/10.17504/protocols.io.bcmriu56>**

**Manuscript citation:**

Chen C, Hao L, Wei W, Li F, Song L, Zhang X, Dai J, Jie Z, Li J, Song X, Wang Z, Zhang Z, Zeng L, Du H, Tang H, Zhang T, Yang H, Wang J, Brix S, Kristiansen K, Xu X, Wu R, Jia H, The female urinary microbiota in relation to the reproductive tract microbiota. GigaByte. 2020.

<https://doi.org/10.46471/gigabyte.9>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** February 17, 2020

**Last Modified:** November 25, 2020

**Protocol Integer ID:** 33169

**Keywords:** dna extraction for human microbe sample, dna extraction, human microbe sample, process of total dna extration, total dna extration, dna, extraction

## Abstract

This protocol is used to clarity the process of total DNA extration for human microbe samples.

## Troubleshooting



- 1 200µl samples were mixed with 500 µl TE buffer and 20 µl of 10 mg/ml lysozyme solution.
- 2 Incubate at 37 °C for 90min. Invert the tube several times to mix every 15 minutes. Centrifuge at 13,000 rcf/min for 10 min at 4 °C and discard the supernatant.
- 3 Add 300 µl of 20 mg/ml proteinase K and 5µl of 10% SDS. Vortex to resuspend cells.
- 4 Incubate at 55 °C in a shaking water bath for 120 min to effect complete lysis. Invert the tube several times to mix every 15 minutes.
- 5 Centrifuge at 13,000 rcf/min for 10 min at 4 °C, then transfer the supernatant to a nuclease-free 2 ml microfuge tube.
- 6 Add 700 µl phenol-chloroform-isoamylalcohol (25:24:1) and vortex to mix well, centrifuge at 13,000 rcf/min for 10 min at 4 °C, then transfer the supernatant to a nuclease-free 2 ml microfuge tube.
- 7 Add 700 µl chloroform-isoamylalcohol (24:1) and vortex to mix well. Centrifuge at 13,000 rcf/min for 10 min at 4 °C, then transfer the supernatant to a nuclease-free 1.5 ml microfuge tube.
- 8 Add 1/10 volume of 3M sodium acetate, 2 µl of 5 mg/ml glycogen and 800 µl cold isopropanol. Incubate at -20 °C for the night.
- 9 Centrifuge at 13,000 rcf/min for 10 min at 4 °C, then discard the supernatant.
- 10 Washed with 800 µl of 70% ethanol. Centrifuge at 13,000 rcf/min for 10 min at 4 °C, then discard the supernatant.
- 11 Wash the precipitate with a second 800 µl of 70% ethanol and centrifuge as above. discard the supernatant.
- 12 Let sit for 15 min to dry the precipitate. Add 1×TE buffer to resuspend the precipitate.



- 13 Add 1µl RNase A to samples and invert tube several times to mix. Incubate at 37 °C for 30 minutes.