



Oct 13, 2022

DNA extraction from mouthwash samples using Qiagen DNeasy PowersoilPro Kit

DOI

<https://dx.doi.org/10.17504/protocols.io.eq2ly7dyrlx9/v1>

Ahmed A Shibl¹, Anique Ahmad¹, Tsedenia Denekew¹, Mamon Abd AlBaqi¹, Aashish R Jha¹

¹New York University, Abu Dhabi

Aashish R Jha: jhaar@nyu.edu



Genetic Heritage Group NYU Abu Dhabi

NYU Abu Dhabi

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.eq2ly7dyrlx9/v1>

Protocol Citation: Ahmed A Shibl, Anique Ahmad, Tsedenia Denekew, Mamon Abd AlBaqi, Aashish R Jha 2022. DNA extraction from mouthwash samples using Qiagen DNeasy PowersoilPro Kit. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.eq2ly7dyrlx9/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: September 18, 2022

Last Modified: October 13, 2022

Protocol Integer ID: 70195

Keywords: dna extraction from mouthwash sample, dna from mouthwash sample, dna extraction, mouthwash sample, using qiagen dneasy powersoilpro kit, qiagen dneasy powersoilpro kit this protocol, dna

Abstract

This protocol describes how to extract DNA from mouthwash samples.

Materials

QIAGEN -DNeasy PowerSoil Pro Kit

Ice

Vortex

Omni Bead Ruptor Elite bead beater

Centrifuge

Liquid N₂

Before start

Set centrifuge to 4°C

Keep CD2 solution on ice or at 4°C

Prepare and label collection tubes, microcentrifuge tubes, MB spin columns, and powerbead tubes



- 1 Thaw the sample on ice for  00:30:00 30m
- 1.1 Transfer the desired sample volume  1-2 mL into a 1.5mL or 2mL eppendorf tube
- 2 Centrifuge the transferred sample at maximum speed ( 17000 x g) at 4°C for  00:10:00 10m
- 3 Discard the supernatant carefully without disturbing the pellet
- 3.1 Repeat Step 1.1 if more volume is needed to see a pellet or a higher yield of DNA is required
- 4 Add  800 µL of CD1 and vortex to resuspend pellet
- 4.1 Spin down briefly
- 5 Transfer entire eppendorf content in a PowerBead Pro Tube
- 6 Secure PowerBead Pro tube onto the bead beater and run at maximum speed ( 17000 x g) for  00:05:00 5m
- 7 Centrifuge the PowerBead Pro tube at maximum speed ( 17000 x g) for  00:01:30 1m 30s
- 8 Transfer the supernatant carefully, without disturbing the pellet, into a clean 2mL microcentrifuge tube
- 9 Add  200 µL of CD2 into the 2mL microcentrifuge tube and vortex for  00:00:10 10s



- 10 Centrifuge the 2mL microcentrifuge tube at maximum speed ( 17000 x g) for  00:01:30 1m 30s
- 11 Transfer the supernatant carefully without disturbing the pellet, into a clean 2mL microcentrifuge tube
- 12 Add  600 µL of CD3 into the 2 mL microcentrifuge tube and vortex for  00:00:10 10s
- 13 Load  650 µL of the lysate onto MB spin columns and centrifuge at maximum speed ( 17000 x g) for  00:01:30 1m 30s
- 13.1 Discard flow-through and repeat step 13 to consume all the lysate
- 14 Place the spin column onto a new collection tube, add  500 µL of EA, and centrifuge at maximum speed ( 17000 x g) for  00:01:30 1m 30s
- 15 Discard flow-through and place the spin column back into the collection tube
- 16 Add  500 µL of C5 onto the spin column and centrifuge at maximum speed ( 17000 x g) for  00:01:30 1m 30s
- 17 Discard flow-through and place the spin column back into the collection tube
- 18 Centrifuge at maximum speed ( 17000 x g) for  00:02:00 and place spin column into 1.5mL eppendorf (elution) tube 2m
- 19 Add  50-100 µL of nuclease-free water to the center of the spin column and leave at room temperature for  00:05:00 5m



20 Centrifuge at maximum speed ( 17000 x g) for  00:01:00 , quantify using Qubit, flash freeze, and store at -20/80 °C

1m