



AUG 14, 2020

🌐 DNA Quantification using the Qubit Fluorometer

Julie Haendiges¹, Ruth Timme¹, Padmini Ramachandran¹, Maria Balkey¹

¹US Food and Drug Administration

GenomeTrakr

Tech. support email: genomeTrakr@fda.hhs.gov



Julie Haendiges

US Food and Drug Administration

DISCLAIMER

Please note that this protocol is public domain, which supersedes the CC-BY license default used by protocols.io.

OPEN  ACCESS



DOI:

dx.doi.org/10.17504/protocols.io.bi8dkhs6

Protocol Citation: Julie Haendiges, Ruth Timme, Padmini Ramachandran, Maria Balkey 2020. DNA Quantification using the Qubit Fluorometer.

protocols.io

<https://dx.doi.org/10.17504/protocols.io.bi8dkhs6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: Jul 31, 2020

Last Modified: Nov 10, 2021

PROTOCOL integer ID: 39909

Keywords: DNA Quantitation,
Qubit dsDNA, Whole Genome
Sequencing, GenomeTrakr

ABSTRACT

This procedure outlines the protocol for quantitation of gDNA for subsequent WGS.

This document applies to all laboratory personnel in the Division of Microbiology (DM) as well as laboratories in the GenomeTrakr Network.

Complete in order:

1. DNA Extraction (Manual DNA Extraction *or* Automated DNA Extraction using the Qiacube)

- Step-by-step procedures to obtain high quality DNA from isolates in TSB for whole genome sequencing

2. DNA Quantitation (included SOP)

- Quantitation of extracted DNA using the Qubit Fluorometer

3. Library Preparation for WGS (Library preparation using Illumina DNA Prep *or* Library Preparation using Illumina Nextera XT)

- Library preparation using NexteraXT *or* Illumina DNA Prep (previously Nextera DNA Flex)

4. Sequencing using Illumina MiSeq

5. Data Quality Checks and NCBI Submission

GUIDELINES

Abbreviations:

BR: Broad Range

DM: Division of Microbiology

HS: High Sensitivity

PI: Principal Investigator

ST1: Standard 1

ST2: Standard 2

MATERIALS

MATERIALS

Qubit dsDNA HS Assay Kit **Thermo Fisher Scientific Catalog #Q32851**

Qubit assay tubes **Thermo Fisher Scientific Catalog #Q32856**

Qubit[®] dsDNA BR Assay Kit **Thermo Fisher Catalog #Q32853**

Supplies:

- Pipette tips, sterile, filtered (assorted volumes)
- 5ml round-bottom polystyrene tubes (Fisher Sci Cat# 14-959-2A or equivalent)

Equipment:

- Qubit 3.0 or 4.0 Fluorometer
- Vortex
- Micropipettes, various volumes
- Microcentrifuge

SAFETY WARNINGS

⚠ Take proper precautions and wear appropriate PPE when handling potentially hazardous chemicals. Ensure that chemicals, spent containers, and unused contents are disposed of in accordance with governmental safety standards.

Qubit dsDNA High Sensitivity and Broad Range Reagent Kit: See

ThermoFisher Scientifics SDSs for additional information

- **HS/BR Reagent 200x in DMSO:** GHS Category 4 for flammability, no data is available on the mutagenicity or toxicity of the reagent, it is known to bind nucleic acid so the same safety precautions should be used as when handling other potential mutagens.
- **HS/BR Buffer:** No GHS classification but may cause irritation with susceptible persons.

BEFORE START INSTRUCTIONS

The buffer and reagent should be stored at room temperature while the standards must be stored at 4oC. The reagent is light and humidity sensitive, it should be stored in the dark and kept in the Ziploc bag with the desiccate pack. When stored as directed, kits are stable for 6 months.

Qubit Working Solution and Sample preparation

- 1 Set up the required number of Qubit tubes for standards and samples. The Qubit dsDNA assays require two standards for calibration. Although, calibrations do not have to be performed each time it is recommended to perform prior to quantifying samples with critical input values.
- 2 Label the tube lids. (Do not label the side of the tube as this could interfere with the sample reading), label tubes for standards as S1 and S2.
- 3 Prepare the Qubit working solution by diluting the Reagent in a 1:200 ratio in Buffer. It is recommended to add an additional sample for overage.

	Total Volume per Sample	10 Samples	20 Samples
Reagent	1 μ l	10 μ l	20 μ l
Buffer	199 μ l	1,990 μ l	3,980 μ l

- 4 Add 190 μ l of working solution to each of the standard tubes.
- 5 Add 10 μ l of each standard to the appropriate tube, then mix by vortexing.
- 6 Add 198 μ l of working solution to each sample tube.

- 7 Add 2 μ l of each sample to the appropriate tube, then mix by vortexing. The final volume should be 200 μ l.
- 8 Allow all tubes to incubate at room temperature for 2 minutes.
- 9 Proceed to Reading Standards and Samples for the appropriate instrument.

Reading Standards and Samples with the Qubit 3.0 Fluorometer

- 10 On the home screen of the Qubit, press **dsDNA**, then select **dsDNA: High Sensitivity** or **dsDNA: Broad Range** depending on the kit being used.
- 11 If calibration is being performed, press **Read Standards** to proceed. (If a calibration is not necessary, you can omit the following steps)
 - 11.1 Insert the tube containing Standard #1 into the sample chamber, close the lid, and press **Read Standard**. When reading is complete, remove the tube.
 - 11.2 Insert the tube containing Standard #2 into the sample chamber, close the lid, then press **Read Standard**. When reading is complete, remove the tube.

11.3 The instrument displays the results of the calibration, for further information on interpretation of the results, refer to the Qubit 3.0 Fluorometer User Guide.

12 Press **Read Samples** to proceed

13 In the **Sample Volume** screen, select the volume of sample added to the tube and the output concentration units (i.e. ng/ μ l)

14 Insert the first sample tube into the sample chamber, close the lid, and then press **Read Tube**.

15 The software displays the results of the sample. The top value (in large font) is the concentration of the original sample. The bottom value is the dilution concentration. Record the concentration of the original sample, remove the tube, and repeat readings and results recording for each additional sample.

Reading Standards and Samples with the Qubit 4.0 Fluorometer

16 The Qubit 4.0 has an on-board Reagent calculator that can be used to quickly determine the correct amount of reagent and buffer required to prepare the working solution.

16.1 On the **Home** screen, press **Reagent Calculator**

16.2 Enter the total number of samples and standards that you will be running.

*Optional: Select **Include overage** if you want to include an additional tube in the total calculated volume.*

16.3 Press **Done** to return to the **Home** screen to run your Qubit Assay

17 On the **Home** screen, select **dsDNA**, then **dsDNA: High Sensitivity** or **dsDNA:Broad Range** depending on the kit being used.

18 If calibration is being performed, press **Read Standards** to proceed. (If a calibration is not necessary, you can omit the following steps)

18.1 Insert the tube containing Standard #1 into the sample chamber, close the lid, and press **Read Standard**. When reading is complete, remove the tube.

18.2 Insert the tube containing Standard #2 into the sample chamber, close the lid, then press **Read Standard**. When reading is complete, remove the tube.

18.3 The instrument displays the results of the calibration, for further information on interpretation of the results, refer to the Qubit 4.0 Fluorometer User Guide.

19 Press **Read Samples** to proceed.

- 20 In the **Sample Volume** screen, select the volume of sample added to the tube and the output concentration units (i.e. ng/ μ l)

- 21 Insert the first sample tube into the sample chamber, close the lid, and then press **Read Tube**.

- 22 The software displays the results of the sample. The top value (in large font) is the concentration of the original sample. The bottom value is the dilution concentration. Record the concentration of the original sample, remove the tube and repeat readings and results recording for each additional sample.

Reference Documents

- 23 [Qubit 3.0 Fluorometer User Guide](#)
[Qubit 4.0 Fluorometer User Guide](#)
[Quick Reference Qubit Assays](#)
[Qubit dsDNA HS Assay Guide](#)
[Qubit dsDNA BR Assay Guide](#)