

May 23, 2023

## DNA quantification with PicoGreen



Forked from [DNA quantification with PicoGreen](#)

DOI

[dx.doi.org/10.17504/protocols.io.6qpvrzb3lmkn/v1](https://dx.doi.org/10.17504/protocols.io.6qpvrzb3lmkn/v1)

Alexander Eiler<sup>1</sup>

<sup>1</sup>University of Oslo

AQUA at UiO



Alexander Eiler

Alex private

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.6qpvrzb3lmkn/v1>

**Protocol Citation:** Alexander Eiler 2023. DNA quantification with PicoGreen. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.6qpvrzb3lmkn/v1>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** June 15, 2018

**Last Modified:** May 23, 2023

**Protocol Integer ID:** 13110

**Keywords:** quantification, DNA, PicoGreen, dna quantification with picogreen, dna quantification, selective for dsdna, dsdna, asymmetrical cyanine dye, dna, ssdna, free dye, dye, sensitive than uv absorbance method, picogreen, rna, uv absorbance method

## Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK


The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to [protocols.io](https://www.protocols.io) is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with [protocols.io](https://www.protocols.io), can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

## Abstract

The Quant-iT™ PicoGreen® dsDNA reagent is a proprietary, asymmetrical cyanine dye. Free dye does not fluoresce, but upon binding to dsDNA it exhibits a >1000-fold fluorescence enhancement. PicoGreen is 10,000-fold more sensitive than UV absorbance methods, and highly selective for dsDNA over ssDNA and RNA. For more information see: <https://www.thermofisher.com/order/catalog/product/P7589>

## Materials

MATERIALS

 Quant-iT™ PicoGreen® dsDNA Assay Kit **Life Technologies Catalog #P7589**

## Troubleshooting



## Before start

**Take out PicoGreen (supplied in the PicoGreen kit) from fridge to thaw, before starting.**  
**Protect it from light.**

PicoGreen is solid at 4°C and thaws at room temperature. The reagent is light sensitive and should be kept wrapped in foil both while thawing and in the diluted stage.

## 1 Prepare TE

Calculate the total volume of 1xTE needed:

- 200 µl per sample
- 100 µl per standard
- Total of 3-400 µl for the DNA standard dilution (depending on the assay you use, see step 2)
- Make enough to count for pipetting losses!!!
- E. g. for high range: 20 samples + 2\* 4 standards + DNA standard the amount is  $20 \times 200 + 8 \times 100 + 300 = 5100 \mu\text{l}$  → make 5500 µl

Dilute from 20xTE stock (supplied in the PicoGreen kit).

- Volume of 20xTE = final volume/20;
- Volume of water = final volume – volume of 20\*TE
- E.g. To make 5 ml: mix 250 µl of 20\*TE with 4750 µl of water.

## 2 Prepare DNA standard

Dilute 100 µg/mL **DNA standard** (supplied in the PicoGreen kit) to working dilution.

- For high range prepare 2 ng/µl: mix 6 µl of DNA stock with 294 µl of 1\*TE
- For middle range prepare 0.5 ng/µl: mix 2 µl of DNA stock with 398 µl of 1\*TE
- For high sensitivity prepare 0.05 ng/µl: mix first 2 µl of DNA stock with 18 µl of 1\*TE (working solution), then mix 2 µl of the working solution with 398 µl of 1\*TE

## 3 Prepare PicoGreen

Calculate the total volume of 1:200 diluted PicoGreen reagent needed.

- For each standard and each unknown sample, a volume of 100 µL will be needed = total number of samples (standard + unknowns) x100 µl + some for pipetting losses.
- e.g. For standard series of 8 and 20 samples (=28 total) prepare  $28 \times 100 \mu\text{l} = 2800 \mu\text{l}$ : → make 3000 µl

Dilute PicoGreen reagent with 1\*TE.

- Volume of PicoGreen = final volume/200;
- Volume of TE = final volume – volume of PicoGreen
- e.g. Mix 15 µl of PicoGreen with 2985 µl of 1\*TE

**KEEP diluted PicoGreen IN THE DARK UNTIL USE!!!**

## 4 Prepare plate - add standards

- Use **NUNC 96-well flat bottom black** plates. Cover with aluminium foil inside the lid.



- Mark samples needed to be measured on the lid and also diagram your plate layout in the table as follow (When possible, measure samples at least in duplicates).
- Plan the standard curve in the microtiter plate.

Table 1 High range standard curve: for 2ul samples, test limit is **0 to 100** ng/ul - for higher concentration dilute your samples.

	Plate Well	volume of TE (uL)	volume (uL) of <b>2ug/mL</b> stock DNA	volume of diluted picog dye (uL)	final DNA concentration in assay (ng/mL)	total DNA in well (ng)
	A1 & A2	0.00	100.00	100	1000	200
	B1 & B2	90.00	10.00	100	100	20
	C1 & C2	99.00	1.00	100	10	2
	D1 & D2	100.00	0	100	0	0

Table 2 Medium range standard curve: for 2ul samples, test limit is **0 to 25** ng/ul

	Plate Well	volume of TE (uL)	volume (uL) of <b>0.5ug/mL</b> stock DNA	volume of diluted pg reagent (uL)	final DNA concentration in assay (ng/mL)	total DNA in well (ng)
	A1 & A2	0.00	100.00	100	250	50
	B1 & B2	90.00	10.00	100	25	5
	C1 & C2	99.00	1.00	100	2.5	0.5
	D1 & D2	100.00	0	100	0	0

Table 3 Low range standard curve: for 2ul samples, test limit is **0 to 2.5** ng/ul

	Plate Well	volume of TE (uL)	volume (uL) of <b>0.05ug/mL</b> stock DNA	volume of diluted pg reagent (uL)	final DNA concentration in assay (ng/mL)	total DNA in well (ng)
	A1 & A2	0.00	100.00	100	25	5



	B1 & B2	90.00	10.00	100	2.5	0.5
	C1 & C2	99.00	1.00	100	0.25	0.05
	D1 & D2	100.00	0	100	0	0

## 5 **Prepare plate - add TE**

- Add 1\*TE to each standard well according to table 1-3.
- Add 95-98 µl of 1\*TE to each sample well. Prepare 2 wells for each sample (duplicates). I.e. if you add 2 µl of sample add 98 µl of 1\*TE in the case of 5 µl sample add 95 µl 1\*TE (see point 6).

## 6 **Prepare plate - add DNA samples**

- Add 2-5 µl of sample to each sample well.


## 7 **Prepare plate - add DNA standard**

- Add diluted DNA standard according to Table 1-3.

## 8 **Prepare plate - add PicoGreen**

- Add 100 µl of diluted PicoGreen to each of the wells (both sample and standard wells) and mix by pipetting 5-10 times.

## 9 Incubate in room temperature in dark for at least 5 minutes and measure with the plate reader (ask for help using the plate reader if using it for the first time).

 00:05:00

## 10 To calculate your DNA concentration:

1. Create a calibration curve in excel or similar from the standards (x axis: measured fluorescens, y axis: total DNA in standard wells in ng from corresponding tables in step 4).
2. Use the equation of the fitted trendline (excel) to calculate the amount of DNA (ng) in your samples.
3. To calculate the DNA concentration of your samples (ng/µl), divide the calculated amount of DNA (ng) by the volume of DNA (µl) that you added in step 6 (2-5 µl).