

Oct 25, 2020

## DNT Detection In Soil

DOI

[dx.doi.org/10.17504/protocols.io.bnxtmfnn](https://dx.doi.org/10.17504/protocols.io.bnxtmfnn)

Zhujun Wei<sup>1</sup>

<sup>1</sup>2020 iGEM NEFU China

2020 iGEM NEFU China



Zhujun Wei

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bnxtmfnn>

**Protocol Citation:** Zhujun Wei 2020. DNT Detection In Soil. protocols.io <https://dx.doi.org/10.17504/protocols.io.bnxtmfnn>

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

We use this protocol and it's working



**Created:** October 25, 2020

**Last Modified:** October 25, 2020

**Protocol Integer ID:** 43731

**Keywords:** dnt detection, soil

## Troubleshooting



- 1 The 1% bacteria (PYB1s-yqjF4-luxCDABE) were cultured in LB containing 0.1% streptomycin (50 mg/mL) for 24 h with aeration at 37°C to reach a final density of  $1 \times 10^8$  cells/mL.
- 2 Preparation of 2% Alginate solution, 1% calcium-chloride solution, experimental soil:dissolve 2g Alginate in 100 mL distilled water (2% w/w) and 1g calciumchloride in 100 mL distilled water (1% w/w). The experiment soil was filtrated prior to use.
- 3 Sterilize the 2% Alginate solution, 1% calcium-chloride solution, (121°C, 15min), and experimental soil(121°C, 1.5h).
- 4 Add 0, 0.25, 0.5and 1mg of DNT powder(0, 250, 500 and 1000mg/kg) to the wells of a 12-well plate.Take 1g of experimental soil and drop it ontothe DNT in the 12-well plate on a clean bench.
- 5 The engineered bacteria were added into 2% Alginate solution at a 1:9 (v/v) ratio.
- 6 Drop 200  $\mu$ L of 1% (w/w) calcium-chloride solution into the 1mL alginate-bacterial suspension droplets (the volume ratio between the Alginate mixture and calcium chloride solution was 5:1). Leave the mix for 1 minute.
- 7 The mixed Alginate gel was tiled on the soil.