

Oct 26, 2020

DNT Detection In Soil

DOI

<https://dx.doi.org/10.17504/protocols.io.bnydmfs6>

Zhujun Wei¹

¹2020 iGEM NEFU China

2020 iGEM NEFU China



Zhujun Wei

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bnydmfs6>

Protocol Citation: Zhujun Wei 2020. DNT Detection In Soil. protocols.io <https://dx.doi.org/10.17504/protocols.io.bnydmfs6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 26, 2020

Last Modified: October 26, 2020

Protocol Integer ID: 43749

Keywords: dnt detection, soil

- 1 The 1% bacteria (PYB1s-yqjF3rd-luxCDABE) were cultured in LB containing 0.1% streptomycin (50 mg/mL) for 24 h with aeration at 37°C to reach a final density of 1×10^8 cells/mL.
- 2 Preparation of 2% Alginate solution, 1% calcium-chloride solution, experimental soil:dissolve 2 g Alginate in 100 mL distilled water (2% w/w) and 1g calciumchloride in 100 mL distilled water (1% w/w). The experiment soil was filtrated prior to use.
- 3 Sterilize the 2% Alginate solution, 1% calcium-chloride solution, (121°C, 15min), and experimental soil (121°C, 1.5h).
- 4 Add 0, 0.25, 0.5 and 1 mg of DNT powder (0, 250, 500 and 1000 mg/kg) to the wells of a 12-well plate.Take 1g of experimental soil and drop it onto the DNT in the 12-well plate on a clean bench.
- 5 The engineered bacteria were added into 2% Alginate solution at a 1:9 (v/v) ratio.
- 6 Drop 200 μ L of 1% (w/w) calcium-chloride solution into the 1 mL alginate-bacterial suspension droplets (the volume ratio between the Alginate mixture and calcium chloride solution was 5:1). Leave the mix for 1 minute.
- 7 The mixed Alginate gel was tiled on the soil.