

Sep 20, 2019

Dot mutation

DOI

<https://dx.doi.org/10.17504/protocols.io.7iehkbe>

宏亮¹

¹Northeast Forest University

2019 iGEM NEFU_China

Tech. support email: shengyiyanwork@gmail.com



宏亮 董

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.7iehkbe>

Protocol Citation: 宏亮 2019. Dot mutation. protocols.io <https://dx.doi.org/10.17504/protocols.io.7iehkbe>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 20, 2019

Last Modified: September 20, 2019

Protocol Integer ID: 27942

Keywords: Mutation, Dot mutation

Safety warnings

 Please wear gloves for the experiment.



- 1 The entire plasmid was amplified reversely by PCR using primers with the fragment sequence that you want to replace.
- 2 The temple plasmids in the PCR process were digested by DpnI enzyme.
- 3 The digested product was transferred into DH5 α . Overnight culture them at 37°C.
- 4 To determine whether the vector was constructed successfully, colony PCR and enzyme digestion were done.
- 5 Select the positive results for sequencing and the final results were obtained.