Oct 24, 2020

## Ot mutation

DOI

dx.doi.org/10.17504/protocols.io.bnwxmffn

Zhujun Wei<sup>1</sup> <sup>1</sup>2020 iGEM NEFU China

2020 iGEM NEFU China

Zhujun Wei





## DOI: dx.doi.org/10.17504/protocols.io.bnwxmffn

Protocol Citation: Zhujun Wei 2020. Dot mutation. protocols.io https://dx.doi.org/10.17504/protocols.io.bnwxmffn

**License:** This is an open access protocol distributed under the terms of the **<u>Creative Commons Attribution License</u>**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working We use this protocol and it's working

Created: October 24, 2020

Last Modified: October 24, 2020

Protocol Integer ID: 43703

## Safety warnings

Please wear gloves for the experiment.

- 1 The entire plasmid was amplified reversely by PCR using primers with the fragment sequence that you want to replace.
- 2 The temple plasmids in the PCR process were digested by DpnI enzyme.
- 3 The digested product was transferred into DH5 $\alpha$ . Overnight culture them at 37°C.
- 4 To determine whether the vector was constructed successfully, colony PCR and enzyme digestion were done.
- 5 Select the positive results for sequencing and the final results were obtained.