

Sep 25, 2022

dsRNAs treatment with RNase T1 and DNase I

DOI

<https://dx.doi.org/10.17504/protocols.io.36wgqj9kyvk5/v1>



Vahid Jalali Javaran¹

¹Département de Biologie, Centre SÈVE, Université de Sherbrooke, Sherbrooke, QC J1K 2R1, Canada



Vahid Jalali Javaran

University of Sherbrooke

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.36wgqj9kyvk5/v1>

Protocol Citation: Vahid Jalali Javaran 2022. dsRNAs treatment with RNase T1 and DNase I. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.36wgqj9kyvk5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: September 16, 2022

Last Modified: September 25, 2022

Protocol Integer ID: 70142

Keywords: dsrnas treatment, samples with rnase t1, rnase t1, total concentration of rna, rna, stranded rna, rnas in the absence, nanopore, treating sample

Abstract

For dsRNA sequencing by nanopore sequencing, this protocol was used. Before treating samples with RNase T1, you should measure the total concentration of RNAs in the samples by using a nanodrop or Qubit device, as RNase T1 has the ability to partially digest double-stranded RNAs in the absence of single-stranded RNA.

Materials



00:00:00

1. DNase I (RNase-free)
2. DNase I Reaction Buffer (10X)
3. RNase T1



Digestion

30m

- 1 Add 10X DNase Buffer with MgCl₂ (final concentration should be 1X).
- 2 Add 50 units RNase T1 per 1µg of total RNA and 1 unit DNase I per 2µg of total RNA
- 3 Incubate at 37 degrees C for 20 min.

Inactivation of enzymes

- 4 cleanup with phenol/chloroform extraction.