

Dec 06, 2019

Version 1

Dye-terminator DNA sequencing V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.97jh9kn>

Diep R. Ganguly¹

¹University of Pennsylvania



Diep R R Ganguly

The Australian National University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.97jh9kn>

Protocol Citation: Diep R. Ganguly 2019. Dye-terminator DNA sequencing. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.97jh9kn>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 06, 2019

Last Modified: December 06, 2019

Protocol Integer ID: 30667

Keywords: Sanger sequencing, Dye-terminator sequencing, sanger sequencing, reactions for sanger sequencing, plasmid dna on the 3130x, genetic analyser, plasmid dna, sequencing reaction, terminator cycling, performing terminator cycling, amplified pcr product, applied biosystem

Abstract

This protocol (based on the BigDye® Terminator v3.1 Cycle Sequencing Kit) is for performing terminator cycling sequencing reactions for Sanger sequencing of amplified PCR products or plasmid DNA on the 3130X genetic analyser (Applied Biosystems).

Attachments



BigDye Terminator v3...

350KB

Materials

MATERIALS

- ⊗ Antarctic Phosphatase - 1,000 units **New England Biolabs Catalog #M0289S**
- ⊗ 96 well PCR Plate Non-skirted **Phenix Research Catalog #MPS-499**
- ⊗ Nuclease-free water (e.g. MilliQ or HPLC grade water)
- ⊗ primers
- ⊗ EDTA
- ⊗ 10 mM dNTPs **Life Technologies Catalog #10297-018**
- ⊗ Ethanol **Merck Millipore (EMD Millipore) Catalog #100983**
- ⊗ BigDye® Terminator v3.1 Cycle Sequencing Kit **Thermo Fisher Catalog #4337454**
- ⊗ Exonuclease I (E. coli) **NEB Catalog #M0293S**
- ⊗ Hi-Di™ Formamide **Thermo Fisher Scientific Catalog #4311320**

Before start

Optimize PCR cycling (if sequencing amplified PCR products) to ensure your reaction produces a single product. Perform gel excision or PCR clean-up with the potential inclusion of incubating with Antarctic phosphatase and Exonuclease 1 to dephosphorylate and degrade unincorporated dNTPs in PCR reactions to prepare templates for DNA sequencing.



Terminator cycling reaction

- 1 Perform sequencing reaction with BigDye Terminator cycling kit and either forward or reverse primers, or both.

Component	Volume (μl)
2.5X Reaction Ready Mix	4
5X BigDye Sequencing buffer	2
20 μM F/R Primer	1
Template (plasmid or cleaned PCR product)	150-300 ng dsDNA or approx 10ng PCR product (see BigDye manual)
Nuclease-free water	to 20 μl

BigDye Terminator Cycling reaction



- 2 Set the following protocol and allow thermal cycler to reach 96 °C:
 1. 1 min at 96 °C
 2. 30 cycles consisting of: 96 °C for 10 seconds, 50 °C for 5 seconds, and 60 °C for 4 min.
 3. Hold at 10 °C.

Purify products

- 3 To the PCR reaction, add 60 µL 100% ethanol and 5 µl 125 mM EDTA.
- 4 Incubate at room temperature for 15 minutes.
- 5 Centrifuge at 4 °C at max speed for 10 minutes.
- 6 Remove as much supernatant as possible, then allow to air-dry in the dark for 10-15 minutes.

Resuspend products and submit for sequencing

- 7 Resuspend the pellet (likely transparent) in 20 µL HiDi Formamide (add 20 µL to any empty wells). Spin down briefly.
- 8 Submit for sequencing on 3130X genetic analyser (Applied Biosystems).