

Jul 12, 2020

Version 5

Dye-terminator DNA sequencing V.5

DOI

<https://dx.doi.org/10.17504/protocols.io.bhxmj7k6>

Diep Ganguly¹

¹University of Pennsylvania

Pogson Group



Diep R R Ganguly

The Australian National University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bhxmj7k6>

Protocol Citation: Diep Ganguly 2020. Dye-terminator DNA sequencing. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.bhxmj7k6>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

Works well for plasmid DNA, optimising sequencing of amplified PCR amplicons.

Created: June 26, 2020

Last Modified: July 12, 2020

Protocol Integer ID: 38605

Keywords: Sanger sequencing, Dye-terminator sequencing, sanger sequencing, reactions for sanger sequencing, plasmid dna on the 3130x, genetic analyser, plasmid dna, sequencing reaction, terminator cycling, performing terminator cycling, amplified pcr product, pcr, applied biosystem

Abstract

This protocol (based on the BigDye® Terminator v3.1 Cycle Sequencing Kit) is for performing terminator cycling sequencing reactions for Sanger sequencing of amplified PCR products or plasmid DNA on the 3130X genetic analyser (Applied Biosystems).

Attachments



[BigDye Terminator v3...](#)

350KB



[wizard-sv-gel-and-pc...](#)

426KB

Materials

MATERIALS

✕ XRN-1 - 100 units **New England Biolabs Catalog #M0338L**

✕ Antarctic Phosphatase - 1,000 units **New England Biolabs Catalog #M0289S**

✕ 96 well PCR Plate Non-skirted **Phenix Research Catalog #MPS-499**

✕ Wizard SV Gel and PCR Clean-Up System **Promega Catalog #A9281**

✕ Nuclease-free water (e.g. MilliQ or HPLC grade water)

✕ primers

✕ EDTA

✕ 10 mM dNTPs **Life Technologies Catalog #10297-018**

✕ Ethanol **Merck Millipore (EMD Millipore) Catalog #100983**

✕ BigDye® Terminator v3.1 Cycle Sequencing Kit **Thermo Fisher Catalog #4337454**

✕ Exonuclease I (E. coli) **NEB Catalog #M0293S**

✕ Hi-Di™ Formamide **Thermo Fisher Scientific Catalog #4311320**



Before start

Optimize PCR cycling (if sequencing amplified PCR products) to ensure your reaction produces a single product. If needed, perform gel excision and clean-up to purify the target DNA fragment. Incubate with Antarctic phosphatase (SAP, AP, or CIP) and Exonuclease 1 to dephosphorylate and degrade unincorporated dNTPs prior to incorporating fluorescent nucleotides in the sequencing PCR (BigDye reaction).



Enzymatic PCR clean-up

1h

- 1 If sequencing a PCR amplified DNA fragment, gel purify target DNA band based on expected fragment size (if multiple bands present). Perform gel purification with Wizard SV Gel and PCR Clean-Up System (Promega, as per attached Manufacturer's instructions) followed by enzymatic clean-up (hydrolyze excess primers and nucleotides) with the following reaction:

Component	Volume (μl)
10X Antarctic phosphatase reaction buffer	1
Antarctic phosphatase	0.5
XRN-1	0.5
Purified DNA fragment	50-150 ng DNA
Nuclease-free water	to 10 μl

Enzymatic clean-up of PCR products

Incubate the above in a thermal cycle for:

1. 37 °C for 30 minutes
2. 80 °C for 15 minutes.

Terminator cycling reaction

- 2 Perform sequencing PCR in PCR tubes (or 96-well plate) with BigDye Terminator cycling kit and forward or reverse primers.

30m

Component	Volume (μl)
v3.1 Ready reaction mix	1
5X Sequencing buffer	1.5
20 μM F/R Primer	0.5



Template (plasmid or cleaned PCR product)	50-150 ng DNA
Nuclease-free water	to 10 μ l

BigDye Terminator Cycling reaction

5x reaction buffer=400 mM TRIS, 10 mM $MgCl_2$

3 Run the following thermal cycling protocol:

1. 1 min at 96 °C
2. 30-40 cycles: 96 °C for 10 seconds, 50 °C for 5 seconds, and 60 °C for 4 min.
3. Hold at 4-12 °C.

4h

Purification

1h 30m

- 4 Transfer PCR reaction to nuclease-free eppendorf tube. To the reaction, add 2.5 μ L of 125 mM EDTA (make sure it touches bottom of tube).
- 5 Add 30 μ l of 100% ethanol, *mix well* (inversion).
- 6 Incubate at room temperature for 15 minutes.
- 7 Centrifuge at 4 °C at max speed for 30 minutes.
- 8 Discard supernatant and add 50 μ l of ice-cold 70% ethanol.
- 9 Centrifuge at 4 °C at max speed for 5 minutes.
- 10 Discard supernatant and allow to air-dry in the dark for >15 minutes.

Prepare for sequencing



- 11 Resuspend the pellet (likely transparent) in 7.5 μ L HiDi Formamide (add to any empty wells). Incubate at RT for 5 minutes then transfer to plate. Spin down briefly.
- 12 Incubate plate at 95 °C for 3 minutes (denature) then place immediately on ice. Spin down briefly.
- 13 Submit for sequencing on 3130X genetic analyser (Applied Biosystems). Keep samples on ice.