

May 07, 2025

Dynabeads Wastewater Virus Enrichment & Total Nucleic Acid Extraction

DOI

<https://dx.doi.org/10.17504/protocols.io.n2bvj9j5nlk5/v1>

Victor Mabasa¹, Natasha Singh¹, Emmanuel Phalane¹, Mokgaetji Macheke¹, Thabo Mangena¹,
Lethabo Monametsi¹, Lebohang Rabotapi¹, Sipho Gwala^{1,2}, Nkosenhle Ndlovu¹, Fiona Els^{1,3},
Sibonginkosi Maposa¹, Said Rachida¹, Kerrigan McCarthy^{1,2}, Mukhlid Yousif^{1,2}

¹National Institute for Communicable Diseases; ²University of Witwatersrand; ³Gauteng City-Region Observatory

Wastewater Genomics S...



Victor Mabasa

National Institute for Communicable Diseases

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n2bvj9j5nlk5/v1>

Protocol Citation: Victor Mabasa, Natasha Singh, Emmanuel Phalane, Mokgaetji Macheke, Thabo Mangena, Lethabo Monametsi, Lebohang Rabotapi, Sipho Gwala, Nkosenhle Ndlovu, Fiona Els, Sibonginkosi Maposa, Said Rachida, Kerrigan McCarthy, Mukhlid Yousif 2025. Dynabeads Wastewater Virus Enrichment & Total Nucleic Acid Extraction. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.n2bvj9j5nlk5/v1>



License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 28, 2025

Last Modified: May 07, 2025

Protocol Integer ID: 119215

Keywords: Wastewater, enrichment, dynabeads , virus enrichment , dynabeads wastewater virus enrichment, total nucleic acid extraction, virus enrichment, kingfisher flex purification system, purified nucleic acid, wastewater sample, wastewater, acid extraction, extraction, digital pcr, extraction process, pcr, virus, generation sequencing

Disclaimer

This protocol was amended from Pub. No. MAN0025695 by appliedbiosystems

Abstract

The Applied Biosystems MagMAX Wastewater Ultra Nucleic Acid Isolation Kit (Cat. No. A52610) is specifically designed for virus enrichment and total nucleic acid extraction from 10 mL wastewater samples. The purified nucleic acids suit various downstream applications, including real-time PCR, digital PCR, and next-generation sequencing. Both the enrichment and extraction processes are automated on the KingFisher Flex Purification System with a 24-deep-well head, using in-house adapted programmes from MagMAX_Wastewater_10mL_Flex24.

Guidelines

General

- Perform all steps at room temperature (20–30°C), unless otherwise noted.
- Clean the work surfaces with RNaseZap to remove nucleases, then wipe the surfaces with 70% to 100% molecular biology grade ethanol to remove additional contaminants.
- Precipitates can form in the Lysis Buffer, Binding Solution, and Wash Buffer if stored below 20°C. If this occurs, warm the reagents at 37°C, then gently mix to dissolve the precipitates. Avoid creating bubbles.

Binding Bead Mix

- Vortex Binding Beads thoroughly before each use.
- Ensure that the beads stay fully mixed within the solution during pipetting.
- Avoid creating bubbles during mixing and aliquoting.
- The binding Bead Mix is very dense so pipet carefully to ensure that the correct volume is added to the sample

Materials

Reagents

Nuclease-free water

Ethanol, 100% (molecular biology grade)

MagMAX Wastewater Ultra Nucleic Acid Isolation Kit (A52610)

Consumables

RNase-Free Microfuge Tubes, 1.5 mL

KingFisher Flex 24 Deep-Well Plate

MicroAmp Clear Adhesive Film

Conical Tubes (50 mL)

Equipment

KingFisher Flex Purification System with 24 deep-well head

Standard laboratory vortex

Pipettes

Automation Scripts

Software

Wastewater_10mL_Virus_Isolation

NAME

Software

Wastewater_10mL_Virus_Extraction

NAME



Before start

- Prepare 80% ethanol using 100% absolute ethanol and nuclease-free water, ensuring a minimum volume of 2 mL per sample.
- Vortex the Binding beads gently to ensure that the beads are fully resuspended.
- Prepare Binding Bead Mix — Combine 500 μ L of Binding Solution with 20 μ L of Binding Beads per sample, preparing enough for the required number of samples plus an additional 10% overage.
- Mix well by inversion, then store at room temperature.



Automated enrichment process

- 1 Set up and label your plates according to the following table:

	Plate Name	Plate type	Contents	Vol per well
	Sample plate 1	24 deep-well	Clarified supernatant	5 000 µL
			Dynabeads	100 µL
	Sample plate 2	24 deep-well	Clarified supernatant	5 000 µL
	Lysis buffer	24 deep-well	Lysis buffer	500 µL
	Tip comb	24 deep-well tip comb	Not applicable	

- 2 Select the appropriate program on the instrument (Wastewater_10mL_Virus_Isolation)
- 3 Start the run, then load the prepared sample and processing plates into position when prompted by the instrument
- 4 While the instrument is running, set up the nucleic acid extraction plates

Automated total nucleic acid extraction process

- 5 Set up and label the Wash, Elution and Tip Comb plates outside of the instrument according to the following table:

6

	Plate name	Plate type	Component	Vol per well
	Wash Buffer 1	24 deep-well	Wash Buffer	1 000 µL

	Plate name	Plate type	Component	Vol per well
	Wash Buffer 2	24 deep-well	80% Ethanol	1 000 µL
	Elution	24 deep-well	Elution Solution	100 µL
	Tip Comb	24 deep-well tip comb		

- 7 When the enrichment process is complete (approximately 20 minutes after starting the run), remove the lysis plate from the instrument and store it safely for nucleic acid extraction. Discard the remaining plates from the instrument
- 8 Add 40 µL of Proteinase K to each enriched sample in the lysis plate
- 9 Invert the Binding Bead Mix tube several times to fully resuspend the beads, then add 520 µL to each sample in the lysis plate
- 10 Load the lysis plate along with Wash 1, Wash 2, Elution, and a new tip comb. Start the run on the instrument and select the appropriate programme (Wastewater_10mL_Virus_Extraction)
- 11 Immediately remove the Elution plate from the instrument at the end of the run and cover it. Alternatively, transfer the eluate to a new tube or plate for final storage. The isolated nucleic acid is ready for immediate use. For storage, keep it at –20°C for up to 6 months or at –80°C for longer than 6 months.

Manual enrichment process

- 12 Follow the manufacturer's protocol available here: https://assets.thermofisher.com/TFS-Assets/LSG/manuals/MAN0025695_MM_Wastewater_Ultra_NA_Isol_Kit_10mL_UG.pdf

Protocol references

This protocol is derived from Pub. No. MAN0025695 by Appliedbiosystems

Acknowledgements

This protocol is derived from Pub. No. MAN0025695 by Appliedbiosystems