

Feb 06, 2019

E. coli Plating Quantification

DOI

dx.doi.org/10.17504/protocols.io.xqhfmt6



Kenneth Schackart¹, Kattika Kaarj¹

¹University of Arizona

481b Laboratory



Kenneth Schackart

University of Arizona

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.xqhfmt6>

Protocol Citation: Kenneth Schackart, Kattika Kaarj 2019. *E. coli* Plating Quantification. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.xqhfmt6>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 03, 2019

Last Modified: February 06, 2019

Protocol Integer ID: 19945

Keywords: plating quantification, plating quantification this protocol, quantification

Abstract

This protocol describes how to serially dilute and plate cultured *E. coli* K-12 in suspension for quantification.

Guidelines

Labcoat and gloves must be worn at all times.


Materials

- Gloves
- Stock *E. coli* K-12 culture
- Pipette and tips
- 2 mL centrifuge tubes
- DI water
- Agar plate
- Inoculation loop
- Parafilm

Troubleshooting

Prepare Serial Dilutions

- 1 Transfer 100 μL of stock solution to 2 mL centrifuge tube.
- 2 Add 900 μL DI water to the same centrifuge tube to make a dilution that is one-tenth the concentration of the stock solution.
- 3 Using 100 μL of the dilution you just made, make another dilution that is one-tenth the concentration of the second solution.
- 4 Repeat 6 times until you have a solution that is diluted by 10^6 .

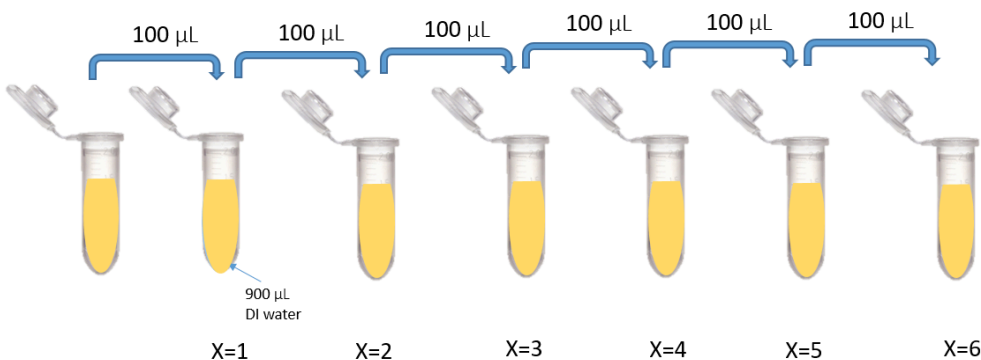

How to perform serial dilutions. X represents the number of 10-fold dilutions performed at each point.

Plate

- 5 Choose a dilution that you feel is appropriate and dispense 100 μL of that dilution on agar plate.
- 6 Evenly spread the cell suspension over the agarose gel with inoculation loop.



Note

This may be different from how you have streaked before. We are wanting the solution evenly spread so that we can count how many colonies are formed after incubation.

- 7 Wrap the agar plate's edge in parafilm.

Note

If you have not worked with Parafilm, ask a TA for help in wrapping your plate.

- 8 Wait 2 minutes, then flip the plate upside-down for incubation.

- 9 Label the plate with:
- Group number
 - Date
 - Dilution used (10^x)

Incubate

- 10 Incubate at  37 °C over night.

Note

Performed by T.A.

Incubate

- 11 Take image of incubated plate.

Note

Performed by T.A.

Include this image in the lab report results section.



Quantify Concentration

- 12 Using the image of your plate, count the number of colonies formed. This can be done by hand or using an image processing software such as ImageJ.
- 13 Calculate the original *E. Coli* concentration using the following equation:

$$C_0 \left[\frac{CFU}{mL} \right] = \frac{\text{colonies counted [CFU]}}{\text{volume dispensed on plate [mL]}} \times \frac{10^x}{1}$$

x represents the number of tenfold dilutions

Include the result of this calculation in the lab report.